

FIELD SERVICE BULLETIN

Protocol 3 Automatic Start

Purpose: Instructions how to use the Protocol 3 control to start a profile at a set time and day.

Accessing the Configuration Menu:

To access the Configuration Menu, and all of its sub-menus, the control must be in its cycle complete or stopped state (not running), if it is not in that state press the button to stop the control before proceeding.

- 1. Press and together from the Select a Mode screen. This should access the Main Menu.
- 2. From the **Main** Menu, use to scroll to **Profile Setup** Menu. Press when menu is highlighted.
- 3. A prompt will appear to enter an unlock code. Enter the correct code with and (the default code is 0010), then press . This should display the **General Configuration** parameter.
- 4. Press when **General Configuration** is highlighted.
 - a. Leaving the **Remote Profile Control** Disabled and Press the to go to **Automatic Start** menu.
- 5. Use the to scroll down to the **Enabled** and Press to enter it.

Entering a New Profile:

- 1. Use the to scroll to down to the **Create a Profile** and Press
- 2. Enter the profile name by changing each letter by using the and and pressing to move across until the end.

MINNEAPOLIS • SHANGHAI • BERLIN • SINGAPORE • HSINCHU



FIELD SERVICE BULLETIN

3.	The Profile Start Trigger menu will be displayed. Use the to scroll to down to the Day and
	Time and Press
	The Profile Start Time will be displayed. using the and change the time and press the to enter it.

- 5. The choice of Days are as follows:
 - a. Monday, Tuesday, Wednesday, Thursday, Friday, Saturday, Sunday, Mon-Fri, Mon-Sat, Sat-Sun, All Week.
- 6. Press when day is highlighted and continue to enter the rest of the profile.

Note: This profile must be started prior to the selected time and day for it to start correctly.