THRIVING ON INNOVATION



LND SERIES OVEN WITH PROTOCOL PLUS™ INSTRUCTION MANUAL

C-187 PN 143377 REVISION I 2/2009

SERVICE AND TECHNICAL SUPPORT

service parts: 1-800-473-7373

international service/main: 1-952-469-8230

service fax: 1-952-469-8193

service@despatch.com

GLOBAL HEADQUARTERS

phone: 1-888-DESPATCH (1-888-337-7282) international/main: 1-952-469-5424

fax: 1-952-469-4513

sales@despatch.com www.despatch.com 8860 207th Street West Minneapolis, MN 55044 USA



Notice

Users of this equipment must comply with operating procedures and training of operation personnel as required by the Occupational Safety and Health Act (OSHA) of 1970, Section 6 and relevant safety standards, as well as other safety rules and regulations of state and local governments. Refer to the relevant safety standards in OSHA and National Fire Protection Association (NFPA), section 86 of 1990.

Caution

Setup and maintenance of the equipment should be performed by qualified personnel who are experienced in handling all facets of this type of system. Improper setup and operation of this equipment could cause an explosion that may result in equipment damage, personal injury or possible death.

Thank you for choosing Despatch Industries. We appreciate the opportunity to work with you and to meet your heat processing needs. We believe that you have selected the finest equipment available in the heat processing industry.

At Despatch, our service does not end after the purchase and delivery of our equipment. For this reason we have created the Service Products Division within Despatch. The Service Products Division features our Response Center for customer service. The Response Center will direct and track your service call to ensure satisfaction.

Whenever you need service or replacement parts, contact the Response Center at 1-800-473-7373: FAX 1-952-469-8193.

Thank you for choosing Despatch.

Sincerely,

Despatch Industries



Standard Products Product Warranty

Products Covered by this Warranty

This warranty (the "warranty") applies to the following Despatch products: LEB, LBB, LAC, LCC, LCD, LLD, RAD, RFD, LND, RTFO, TAD, TFD, PR, PN, PW, PTC and the following Ransco products: RTH, RTS, 900 Series.

Parts and Materials

Despatch warrants all parts and materials to be free from defects in material and workmanship for a period of:

- Five (5) years from date of shipment for laboratory oven electric heaters.
- Three (3) years from the date of shipment for Protocol Plus and DES 2000 temperature controllers; and
- One (1) year from the date of shipment, or 2,000 hours of operation, whichever occurs first, for all other components of products covered by this Warranty.

During the applicable Warranty period, Despatch will repair or replace, at Despatch's option, parts and materials covered by this Warranty.

Labor

During the first 90 days of the Warranty period, Despatch will pay labor costs incurred to remove defective parts and materials, and to reinstall repaired or replacement parts or materials; provided, however, that Despatch's obligation to pay such labor costs shall be subject to the limitation that the removal and/or reinstallation service must be performed by a Despatch-authorized technician from Despatch's worldwide network of factory-trained professionals at a location within the contiguous United Sates.

Transportation Costs

All transportation costs to transport defective parts or materials to Despatch and to transport repaired or replacement parts or materials to Customer shall be the responsibility of Despatch.

Terms and Conditions

This Warranty shall be deemed valid and binding upon Despatch if and only if the Customer:

- Installs, loads, operates, and maintains the covered product supplied hereunder in accordance with the instruction manual provided upon delivery and product labeling affixed to the subject equipment;
- 2. If applicable, follows the Emergency Procedure set forth in this Warranty; and
- Contacts Despatch's Helpline at 1-800-473-7373 for assistance in diagnosing and troubleshooting the problem immediately upon discovering any damage or malfunction

Despatch's reasonable determination as to whether a repair, replacement, or service is covered by this Warranty shall be final and binding.

Exclusions

This Warranty DOES NOT cover,

1. Damage or malfunctions, or expenses incurred in the process of diagnosing and/orrepairing damage or malfunctions, resulting from any of the following: operator

error, misuse, abuse, inadequate preventative maintenance, normal wear and tear, service or modifications by other than Despatch authorized technicians, use of the covered product that is inconsistent with the operation manual or labeling, acts of nature (including, without limitation, floods, fire, earthquake, or acts of war or civil emergency), internal or external corrosion, or non-conforming utilities (including, without limitation, electrical, fuel supply, environmental and intake/exhaust installations);

- Repair or replacement of parts or materials designed and intended to be expendable or consumable; refrigerants, filters, lamps;
- 3. Routine maintenance or
- Labor costs incurred for troubleshooting, diagnostics, or testing (except for testing required to verify that a covered defective part or material has been repaired).

Limitations of Liability

Despatch shall not, in any event, be liable for indirect, special, consequential, incidental, or punitive damages or penalties of any kind, including without limitation loss of revenue, profits or business opportunities resulting from interruption of process or production. In no event shall Despatch be liable for damages in excess of the amounts paid by Customer to Despatch with respect to the applicable product(s). This Warranty does not cover, and Despatch shall not be liable for any losses, costs, damages or expenses resulting from delays in diagnosing or repairing the products, supplying or obtaining replacement parts of materials, strikes, labor stoppages or shortages, fires, accidents, government acts or regulations, or any other causes beyond the control of Despatch.

Non-Compliance By Customer

Despatch reserves the right to suspend and withhold service under this Warranty in the event of non-compliance by the Customer to any terms and conditions of this Warranty or the applicable purchase order or invoice. Further, Despatch shall not be liable for any loss of production, expenses, and inconveniences incurred due to such suspension.

<u>Customer Furnished Equipment Warranty Limitation</u>

This Warranty does not cover diagnosis or repairs of defects in or caused by, lack of performance of, or fitness for purpose of customer-supplied parts or equipment unless specifically noted in the Despatch written order acceptance confirmation.

Performance Commitment

Despatch provides no guarantee of process performance or fitness for purpose, unless specifically noted otherwise in Despatch written order acceptance confirmation. Despatch is providing equipment with design parameters specific only to its equipment.

Procedure Upon Discovery of Defects and Emergencies

In the event Customer becomes aware of any defect in the application products, Customer must immediately: (a) shut off fuel or energy supply (gas and electricity); (b) call for emergency assistance, if needed, and (c) notify Despatch Service.

THE REPRESENTATION AND WARRANTIES SET FORTH HEREIN ARE EXCLUSIVE AND IN LIEU OF, AND CUSTOMER HEREBY WAIVES AND DISCLAIMS RELIANCE UPON, ALL OTHER REPRESENTATIONS AND WARRANTIES OF EVERY KIND WHATSOEVER WHETHER EXPRESS OR IMPLIED, OR ARISING BY OPERATION OF LAW OR IN EQUITY, OR BY COURSE OF PERFORMANCE OR DEALING OR USAGE OF TRADE, INCLUDING, WITHOUT LIMITATION, ANY IMPLIED WARRANTIES OF MERCHANTABILITY OR OF FITNESS FOR A PARTICULAR PURPOSE.

THIS WARRANTY IS PERSONAL TO THE CUSTOMER AND MAY NOT BE TRANSFERRED OR ASSIGNED. ALL LIMITATIONS HEREUNDER, HOWEVER SHALL BE BINDING ON ALL SUCCESSORS AND ASSIGNS OF CUSTOMER.

Despatch Industries Advantage Service Assurance Program (ASAP)

CONTACT: DESPATCH SERVICE AGREEMENTS SPECIALIST at 800-473-7373 or 952-469-8230 or e-mail: service@despatch.com

Despatch continues to deliver exceptional products backed by a strong sense of responsibility and drive for long term customer satisfaction. Your partnership with Despatch can offer even higher value through your subscription to one of Despatch's Advantage Service Assurance Program(ASAP).

Warranty

Despatch's exclusive, comprehensive service programs start with the 1 year parts only warranty which is described on the other side of this document. This warranty can be expanded immediately to meet your most stringent service needs. Despatch Service Products Group will be able to answer your service questions and provide a quotation for the immediate expansion of your product warranty. Call 800-473-7373 or 952-469-8230; or e-mail service@despatch.com.

Immediate Service Response

The key to an effective service program is response. Wherever your location, Despatch is only a phone call away. Our U.S. and Canadian customers can reach Despatch at 1-800-473-7373. Worldwide customers can call 1-952-469-8230 or FAX 1-952-469-8193. Our Customer Service Technicians have over 150 years combined experience and access to detailed design and manufacturing documentation specific to your Despatch unit(s). This exacting level of service is a benefit only Despatch can provide and means that you can expect speedy, accurate and the most cost effective response.

Field Service Network

A worldwide network of factory trained Service Professionals is available to support your Despatch equipment. From routine repair to certified instrument calibration, the Despatch service network is positioned to respond to your needs. As a manufacturer of custom equipment, our service programs are customized to meet your specific needs regarding:

- 1. Service scope
- 2. Response time
- 3. Preventive maintenance frequency and content
- 4. Payment method

Sustained Service Support

At Despatch, long term customer satisfaction means more than just responding quickly and effectively to our customers' service needs. It means offering comprehensive customer support well beyond the scope and duration of our initial warranty. Despatch offers two basic service packages which are customized to each individual customer's need. These service packages are titled Full Service and Preventive Maintenance Plus+ service agreement products. Each is unique in the industry and offer the following benefits:

- 1. Priority response for minimum production interruption
- Preventive maintenance for longer product life
- 3. Discounts on parts and services
- 4. Various payment plans to ease budgeting and recording expenses
- 5. Reduce purchase ordering costs

PREFACE

This manual is your guide to the Despatch oven. It is organized to give you the information you need quickly and easily.

The INTRODUCTION section provides an overview of the Despatch oven.

The THEORY OF OPERATION section details the function and operation of assemblies and subassemblies on the Despatch oven.

NOTE:

Read the entire INTRODUCTION and THEORY OF OPERATION <u>before</u> installing the oven.

The INSTRUCTIONS section provides directions on unpacking, installing, operating and maintaining the Despatch oven.

The APPENDIX section contains Special Instructions for operating the control instrument, a Troubleshooting Table, a list of Accessories and a Warranty.

An efficient way to learn about the oven would be to read the manual while working with the corresponding oven control system. This will give you practical hands-on experience with information in the manual and the oven.

WARNING:

Failure to heed warnings in this instruction manual and on the oven could result in personal injury, property damage or death.

Before operating the equipment, be sure you understand all of the technical information contained in this manual. Information skipped, not understood or misunderstood could create the possibility of operating the equipment in an unsafe manner. This can cause damage to the oven or personnel or reduce the efficiency of the equipment.

Revision B: Corrected Sensor Calibration Page instruction, pages 38-39

Revision C: Modified per Rev C Protocol Plus software

Revision D: Update Despatch Warranty

Revision E: Update to Software Revision 4.0

Revision F: Revised Protocol Plus times. Updated Despatch address

Revision G: Revised warranty

Revision H: Revised Unpacking and Inspection Instructions

Revision I: Updated Control section.

TABLE OF CONTENTS

INTRODUCTION	1
Special Features	1
Specifications	
Dimensions	
Capacities	
Temperature	
Power	
THEORY OF OPERATION	
Inert Atmosphere Theory	6
OVEN INSTRUCTIONS	7
Unpacking and Inspection	7
Set-up	8
Operating	11
Loading the Oven	
Pre-Startup Checklist	
Startup	12
Automatic Nitrogen and Water Cooling	
CONTROL	
Theory of Control Operation	
Operating Modes	
Setup Mode	
Fast Start Mode	
High Limit	
Indicators	
Displays	
Key Functions	
Outputs	
Relay (Continued)	
Communication	
Optional Software	22
INSTRUCTIONS	23
Start-Up	23
Operation	
Manual Mode	
Timer Mode	25
Profile Mode	26
Auto Start Mode	26
Setup Mode	27
Instructions for Setup Mode Pages	28
Program Page	28
Profile #	30
Sample Profile	31
Auto Start Page	
PID Page	
Control Page	
Communication Page (optional)	
Real Time Clock Page	
Relay Outputs Page (optional)	

Test Page	38
Zone Calibration Page	
Sensor Calibration Page	
Enable Page	
Digital Inputs (optional)	
Error Messages and Alarms	
Quick Reference and Default Values	
Technical Specifications	
MAINTENANCE	54
Checklist	54
Tests	
Nitrogen Supply	55
Replacement	
Parts	
Protocol Plus™ Instrument	57
Heater Unit	
Fan Motor	
TROUBLESHOOTING	60
APPENDIX	62
Temperature Scale Conversion (C/F)	
Optional MRC5000 Recorder Setup	

INTRODUCTION

This section provides an overview of the Despatch LND Series forced air oven. The LND Series Ovens have the most effective heat distribution and the fastest processing time of any lab oven their size. Air is discharged from the left side wall of the oven and circulates through the chamber.

Special Features

- The sturdy construction and three inch insulation of the Despatch LND Series ovens contribute to excellent temperature uniformity.
- Other special features include the following:
- Unique Despatch design to combine higher fan volume of forced recirculated air with a system of perforated stainless steel walls for the ultimate in temperature uniformity.
- Welded double wall construction and fiberglass insulation to reduce heat loss.
 Silicone rubber gaskets further minimize heat leakage.
- · Rapid response heater.
- Scratch-resistant baked enamel exterior and stainless steel interior for easy cleaning.
- Purge and maintain valves and flowmeters for creating an inert gas atmosphere in the oven chamber
- · Water cooling coil for dissipating heat from the oven chamber

Specifications

Dimensions

Model	0	nber S n (cm)	Size	Capacity feet ³ Overall Size in (cm)			Maximum Number of		
	W	D	Н	(liters)	W	D	Н	Shelf Positions	
LND 1-42	20 (50)	18 (46)	20 (51)	4.2 (120)	33 (84)	41 (104)	45 (114)	9	
LND 2-11	38 (96)	20 (50)	26 (66)	11.0 (310)	54 (137)	40.5 (103)	63 (160)	6	
LND 2-24	48 (122)	24 (61)	36 (91)	24.0 (680)	66 (166)	44.5 (113)	74 (188)	9	

Capacities

Model		LND 1-42	LND 2-11	LND 2-24
Maximum Load	Lbs	150	180	300
Maximum Shelf Load	Lbs	30	60	100
Recirculating Fan	CFM	300	1,000	1,500
	HP	1/4	1	2
Approximate Weight Net	Lbs.	380	745	1,065
	KG	173	339	484
Approximate Shipping Weight	Lbs.	510	1,045	1,465
	KG	232	475	666

Temperature

Model		LND 1-42	LND 2-11	LND 2-24
Time to Temperature (approximate minutes with no load)	25°C - 100°C 25°C - 200°C 25°C - 343°C	6 25 56	5 14 32	3 12 30
Recovery Time Door Open 1 Min. (approximate minutes with no load)	100°C 200°C 343°C	<1 3 10	<1 2 7	<1 2 4
Temperature Uniformity at	100°C* 200°C* 343°C*	±1°C ±2°C ±3°C	±1°C ±2°C ±3°C	±1°C ±2°C ±3°C
Minimum Operating Temperature** (Approximate w/20EC ambient)	with water cooling without water cooling	55°C 97°C	55°C 100°C	60°C 110°C
Control Stability		±0.3°C	±0.3°C	±0.3°C
Repeatability		±0.5°C	±0.5°C	±0.5°C
Cooldown Time** (approximate minutes with no load)	343 - 60°C	30	35	45

^{*} Figures are based on actual tests in an empty oven. Uniformity can vary slightly depending on unit and operating conditions.

^{** 60°}F Supply water temperature at 120 GPH for LND1-42, 210 GPH for LND2-11 and LND2-24. Minimum operating temperature is also affected by inert gas flow rate and temperature.

Power

Line voltages may vary in some geographical locations. If your line voltage is much lower than the oven voltage rating, warm up time will be longer and motors may overload or run hot. If your line voltage is higher than name plate rating, the motor may run hot and draw excessive amps.

If the line voltage varies more than 10% from the oven voltage rating, some electrical components such as relays, temperature controls, etc. may operate erratically.

Power Requirements

Model	Volts	Amps	Hertz	Phase	Heater KW
LND 1-42	240*	15.6	50/60	1	ω ο
LND 2-11	208	29.8	60	3	
LIND Z II	240	26.1	60	3	9
LND 2-24	480	13.1	60	3	9
	208	52.4	60	3	16
	240	46.0	60	3	16
	480	23.0	60	3	16

^{*} Oven designed for 240 volts (see name plate on oven) will operate satisfactorily on a minimum of 208 volts, but with a 25% reduction in heater power. If your power characteristics are lower, contact Despatch Industries. An option is available to regain the full heater power when operating on 208V.

THEORY OF OPERATION

This section details the function and operation of assemblies and subassemblies on the Despatch LND Series Ovens. These ovens have the most effective heat distribution system and the fastest processing time of any lab ovens their size. They are especially useful for testing, preheating, sterilizing, drying, aging and curing as well as other production applications. Horizontal airflow with precision digital control delivers uniform, fast processing. The overall result is efficient productivity under strenuous conditions.

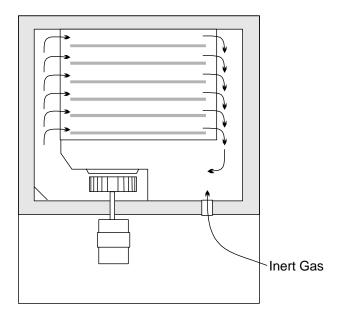


Figure 1. Horizontal Airflow

The unique Despatch computerized design, moves forced convected heat through perforated stainless steel walls. The air is recirculated with a high volume fan. Despatch LND Series Ovens employ higher volume fans than any competitive model. The chamber can be densely loaded without interfering with the process. The LND series air delivery temperature is within 1/2° C of the number appearing on the digital display. Nitrogen intake is regulated by separate purge and maintain panel-mounted flowmeters. The cooling water flow is controlled by an adjustable valve and monitored by a flowmeter located on the rear of the oven.

Inert Atmosphere Theory

The goal of the inert atmosphere oven is to reduce oxygen concentration levels to an acceptable level. This is accomplished by injecting an inert gas into the sealed chamber. As the flow rate of inert gas into the chamber is increased, the chamber becomes pressurized and gas is exhausted out the pressure relief valve located on the back of the oven. Since oxygen is repelled by inert gas flow, the overall oxygen concentration in the chamber is reduced.

However, there is a limit to the desired inert gas flow rate into the chamber. To minimize the oxygen concentration level, you must minimize the number of exhaust outlets. The only desired outlet is through the pressure relief valve. If the inert gas flow rate is too high, other exhaust outlets will be formed (usually the door seal). In this situation oxygen can penetrate back through the leak or outlet even though the chamber is pressurized. This process is called diffusion. As diffusion is minimized, oven performance is maximized.

WARNING:

This oven is not designed for use with flammable material. If your process involves flammable material see the LFD series.

OVEN INSTRUCTIONS

The INSTRUCTIONS section provides directions for unpacking, installation, operation and maintenance of the LND Series oven.

Unpacking and Inspection

Remove all packing materials and thoroughly inspect the oven for damage of any kind that could have occurred during shipment.

- See whether the carton and plastic cover sheet inside carton are still in good condition.
- Look at all outside surfaces and corners of the oven for scratches and dents.
- Check the oven controls and indicators for normal movement, bent shafts, cracks, chips or missing parts such as knobs and lenses.
- Check the door and latch for smooth operation.

If there is damage that may have occurred during shipment, follow these instructions.

- 1. Contact the shipper immediately and file a written damage claim.
- 2. Contact Despatch Industries to report your findings and to order replacement parts for those that were damaged or missing.
- 3. Send a copy of your filed damage claims to Despatch.
- 4. Next, check to make sure you have received all the required materials. Your shipment should include:
 - One (1) Despatch oven,
 - One (1) Instruction manual,
 - One (1) Warranty card,
 - Two (2) Shelves
- 5. If any of these items are missing from the packaged contents, contact Despatch Industries to have the appropriate materials forwarded to you.
- 6. Finally, to protect the warranty on your new LND Series Oven, complete the warranty card and mail it to Despatch within 15 days after receipt of the equipment.

Set-up

1. Select the location for installing your oven.

The oven must have a minimum 2 inch clearance in the rear to provide proper ventilation. The oven may be placed next to another cabinet, or next to another oven, with a 3 inch clearance to insure proper ventilation. Make sure oven is level and plumb; this will assure proper heat distribution and operation of all mechanical components.

WARNING:

Do not use the oven in wet, corrosive or explosive atmospheres unless this oven is specifically designed for a special atmosphere.

- Connect the nitrogen supply line to the inlet located on the oven rear panel and marked nitrogen supply. The nitrogen supply to the oven must not exceed 80 PSI. (40 PSI on LND 1-42)
- 3. Check for leaks in the nitrogen supply line.
 - a. Turn the three-way nitrogen valve switch to the OFF position.
 - b. Turn the PURGE flowmeter knob clockwise to the OFF position.

WARNING:

Use only inert gases.

- c. Turn the MAINTAIN flowmeter knob clockwise to the OFF position.
- d. Open tank valve on the nitrogen supply.
- e. Set pressure regulator to about 40 PSI.
- f. Check the nitrogen plumbing for leaks with a soapy water solution.

As nitrogen gas is odorless, all leaks should be stopped to prevent the possibility of suffocation in a small work area. Gas from a nitrogen leak might leak might displace much of the oxygen in the atmosphere.

4. If cooling is required, install water connection for cooling coils.

WARNING:

All grounding and safety equipment must be in compliance with applicable codes, ordinances and accepted safe practices.

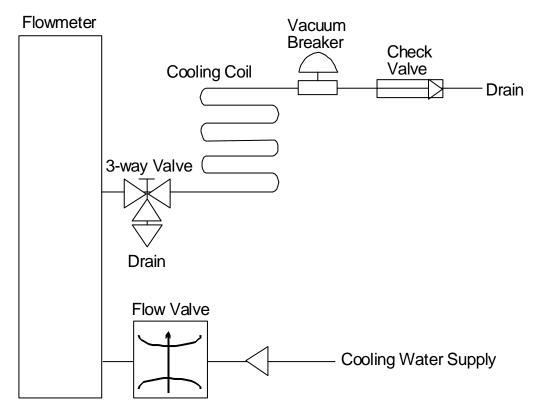


Figure 2. Recommended Water Connection on Despatch LND Series Ovens

- a. Make drain connections at the check valve and 3-way valve.
- b. Connect cooling water supply to the flow valve.
- c. Open 3-way valve and adjust flow control valve for maximum flow.
- d. Check for leaks.

WARNING:

NEVER use the water cooling coil without a gas flow into the chamber. Not using inlet gas during cooldown creates a vacuum pressure which could cause extensive damage to the oven.

CAUTION:

Design the drain system to prevent operator injury from high temperature or pressure buildup. Piping must be able to withstand short periods of up to 650°F (343°C) temperatures. Drain lines should be insulated or warning labels installed that a hazard exists.

WARNING:

Never allow drain to be plugged as a hot oven will generate a small amount of steam when the water is first turned on. Steam burns.

- e. Adjust flowmeter for desired flow rate. Consult the following chart for cooldown rates.
- f. Close 3-way valve. Make sure check valve is working properly by monitoring water flow back through the cooling coil and out the drain port of the 3-way valve.

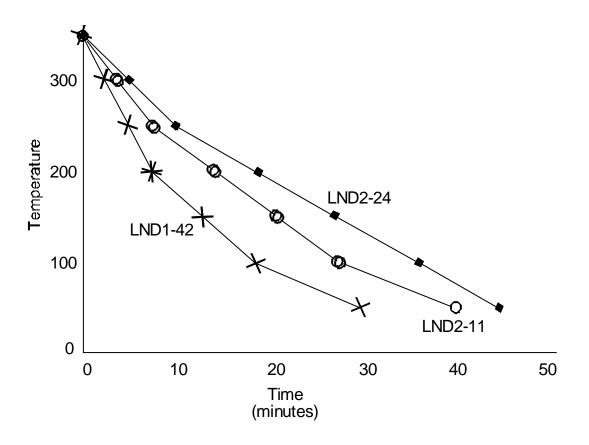


Figure 3. Cool Down Rate with Water Cooling Coil (unit without a load).

NOTE:

Cooldown rates are approximate. Load mass, inert gas flow and exhaust rates affect cool down times.

- 5. Identify correct power source indicated on the specification plate.
- 6. On larger units, remove the control panel and the filter panel to expose the equipment mounting area.
- 7. Hard wire oven directly to the electric supply. On larger units the line connection is located on the equipment mounting panel.

Operating

Users and operators of this oven must comply with operating procedures and training of operating personnel as required by the Occupational Safety and Health Act (OSHA) of 1970, Section 5 and relevant safety standards, and other safety rules and regulations of state and local governments. Refer to the relevant safety standards in OSHA and National Fire Protection Association (NFPA), Section 86 of 1990.

WARNING:

Do not use oven in wet, corrosive or explosive atmospheres unless this oven is specifically designed for a special atmosphere.

Loading the Oven

Despatch Industries cannot be responsible for either the process or process temperature used, or for the quality of the product being processed. It is the responsibility of the purchaser and operator to see that the product undergoing processing in a Despatch oven is adequately protected from damage.

Carefully following the instructions in this manual will help the purchaser and operator in fulfilling that responsibility.

When loading the oven avoid spills of anything onto the heater elements or onto the floor of the oven. Do not place the load on the oven floor plate. This may cause the

load to heat unevenly and the weight may cause shorting out of the heater elements. Use the shelves provided.

The two shelves are designed to be pulled out about half way without tipping. The support capacity of the shelves is listed in the Capacities Table in the Specifications section in this manual. Do not overload the shelves.

WARNING:

Do not use flammable solvent or other flammable material in this oven. Do not process closed containers of any substance because they may explode under heat.

Distribute the workload evenly so that airflow is not restricted. Do not overfill your oven. The workload should not take up more than two-thirds of any dimension of the inside cavity.

Pre-Startup Checklist

- Know the system. Read this manual carefully. Make use of its instructions and explanations. Safe, continuous, satisfactory, trouble-free operation depends primarily on your understanding of the system and your willingness to keep all parts in proper operating condition.
- Check line voltage. Voltage must correspond to nameplate requirements of motors and controls. Refer to the section on power connections in the INTRODUCTION of this manual.
- <u>Fresh air and exhaust.</u> Do not be careless about restrictions in and around the fresh air and exhaust openings and stacks. Under no condition permit them to become so filled with dirt that they appreciably reduce the air quantity. The proper ventilation clearances should be fulfilled at all times. Refer to the Set-up instructions in this manual.

Helpful hints

For drying ovens, open vent to prevent buildup of moisture.

For sample heating, close the vent when no ventilation is required.

Startup

For fastest oven heat-up time, close the fresh-air vent. After the desired temperature is reached, the vent may be adjusted as needed.

- 1. Start Fan.
 - a. Open oven door.
 - b. Press Power switch to the On position. You will hear the recirculating fan start.
 - c. Shut oven door.
 - d. Check that the control display turns on.
- 2. Adjust nitrogen flow rate.
 - a. Determine the desired oxygen level.
 - b. Use the following chart to determine the time to achieve the desired oxygen level.

c. Turn the nitrogen valve switch to the PURGE position for the time to achieve oxygen level as determined in step 2b.

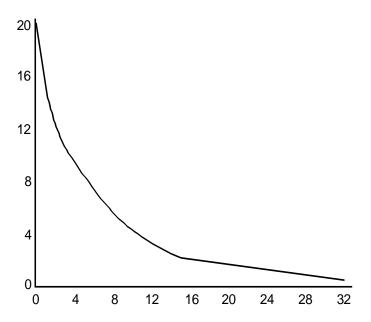


Figure 4. Oxygen Level vs. Time at Maximum SCFH Purge Flow

- d. When the purge is complete, turn the nitrogen valve switch to the MAINTAIN position.
- e. Use the charts on the next page to determine the MAINTAIN flowmeter value to maintain the oxygen stabilization level.
- f. Adjust the MAINTAIN flowmeter to the value determined in step 2e.
- 3. Operate the temperature control as desired by following the control operation instructions that follow.

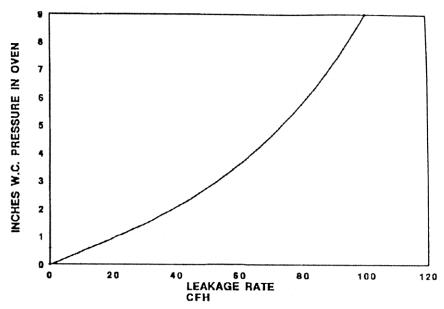


Figure 5. Nitrogen Leakage Rate vs. Oven Pressure

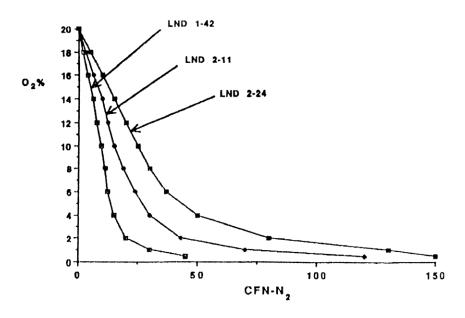


Figure 6. Oxygen Stabilization Level vs. Flow of Nitrogen

Automatic Nitrogen and Water Cooling

Your oven may be fitted with optional automatic water cooling and/or nitrogen control features. These features replace the manual on/off valves with solenoid valves which can be operated either by switches in the control panel or programmed event outputs on the temperature control instrument.

The additional switches in the control panel have AUTO and MANUAL positions. Placing the switch in MANUAL energizes the appropriate solenoid valve. Leaving the switch in the AUTO position allows the solenoid valve to be operated by the control event outputs.

Refer to the following section on the control instrument for more information on programming the event outputs. The standard configuration for these optional features is for event relay 1 to operate water cooling, event relay 3 to operate nitrogen purge, and event relay 4 to operate nitrogen maintain. In addition, if the optional door lock feature is also installed, it will be in the locked position any time the nitrogen maintain solenoid is energized.

Typically, the control is programmed so that the purge and maintain valves are energized at the beginning of a program in a segment of long enough duration to purge the oven (see the oxygen level chart discussed earlier). The next program segment then de-energizes the purge valve while leaving the maintain valve energized. The maintain valve should be left energized for as long as the nitrogen level is desired to be maintained.

The water cooling solenoid valve is energized in the last segment(s) of the program to bring the chamber temperature down to a safe unloading temperature.

CONTROL


The special features of the Protocol PlusTM control include:


- PID tuning
- Ramp/Soak programming of up to 64 segments
- Segment looping and profile linking
- Built-in manual reset high limit control
- Built-in process timer
- Dedicated LED display for process temperature
- Multi-purpose two-line LCD display with backlight
- Auto-tuning
- Security access
- Process temperature retransmission signal
- Digital inputs for remote profile control
- Real time clock
- Optional relay outputs for events, alarms, or end-of-cycle signal
- Optional RS232/RS422/RS485 MODBUS communications

Theory of Control Operation

The Protocol Plus is a modular microprocessor based digital temperature controller. The Protocol Plus operates as a dual functioning controller/high limit instrument. The control portion utilizes a time-proportioning voltage signal to control heating devices with minimal temperature fluctuations.

The high limit portion protects the product and/or the oven from overheating. If the product being processed has a critical high temperature limit, the high limit setpoint should be set to a temperature somewhat below the temperature at which the product could be damaged. If the product does not have a critical high temperature limit, the high limit setpoint should be set 5 to 15 degrees higher than the maximum programmed setpoint at which the oven will operate.

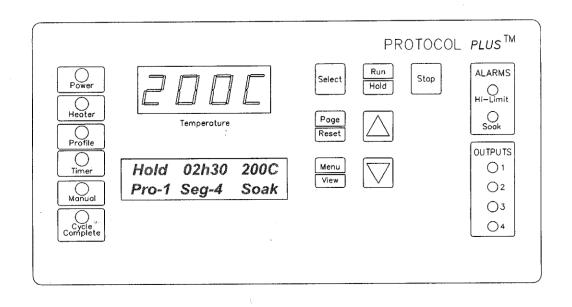

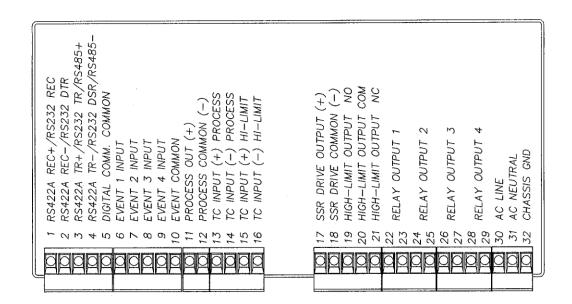

Protocol Plus Faceplate and Wiring Diagram

Operating Modes

The Protocol Plus control has five modes of operation available:

Stopped Mode: All control and relay outputs are off. Stopped Mode is integrated

into each of the following four modes of operation.

Manual Mode: Control operates as a single setpoint control until Stopped mode is

accessed

Timer Mode: Control operates as a single setpoint control until preset time period

has expired.

Profile Mode: Control operates as a ramp/soak profiling control until the end of

the profile. 8 profiles are available with up to 8 ramp/soak segments

in each profile.

Auto Start Mode: Control may automatically start Manual, Timer, or Profile mode

based on a preset time and day.

The optional event outputs can be utilized during Manual, Timer, or Profile modes.

Setup Mode

The control has a Setup Mode which provides access to control configuration and programming of profiles. The Setup Mode contains ten separate electronic Pages where the configuration and programming parameters (Menu items) are found. The Setup Mode Pages are described in detail elsewhere in this manual.

Fast Start Mode

The Protocol Plus control has the ability to automatically start an operating mode when power is applied. This feature may be useful if the same mode of operation is used everyday. The user can turn on the power and the oven will start the desired process automatically. The Fast Start Mode is controlled by the Power-Up Start parameters on the Control page (see Setup Mode).

High Limit

The control has an integrated high limit function which will disable the heater output when tripped. If the high limit does trip, the relay will need to be manually reset. When the high limit relay is tripped, the Hi-Limit indicator will be lit. Allow the oven to cool several degrees (or increase the high limit setpoint) and then press the Reset key. The indicator will turn off.

High-Limit temperature readout is provided on LCD Line #2 in all Modes (Stop, Run, Hold, and Standby) except Setup Mode. High-Limit temperature is displayed for 10 seconds, alternating with current Mode and Status display for 10 seconds.

The control will not allow the high limit setpoint to be set below the current setpoint value.

Indicators

The Protocol Plus control has 12 indicating LEDs that provide operational information to the user.

- Power LED: Indicates that power is supplied to the instrument.
- Heater LED: Indicates that the heater output is active.
- **Profile LED:** Indicates that the Profile Mode is in operation.
- **Timer LED:** Indicates that the Timer Mode is in operation.
- Manual LED: Indicates that the Manual Mode is in operation.
- Cycle Complete LED: Indicates that the control is in Stopped mode.
- Hi-Limit Alarm LED: Indicates that the high limit relay has tripped (de-energized).
- Soak Alarm LED: Indicates that the guaranteed soak deviation is in alarm condition.
- Outputs 1 through 4: Indicate that the optional relay outputs are in the ON state. These outputs may be configured as timed event outputs, process temperature trip point outputs, alarm outputs, or as an end of cycle relay output. The ON state can be configured as energized or de-energized.

Displays

The Protocol Plus control has two displays. A dedicated LED upper display shows the oven temperature. A two-line LCD lower display provides information on control status, high limit temperature, and allows changes to be made to the control settings.

Key Functions


The Protocol Plus control has seven keys that provide operation.

- **Select key:** Press to select mode of operation. In Setup Mode, to select profile number or relay. In Profile/Run Mode, press simultaneously with the UP key to advance a segment.
- Run/Hold key: Press to activate a mode of operation. If a Profile (or Timer)
 Mode is running, pressing the Run/Hold key will place the Profile (or Timer) in
 Hold status. A subsequent press will resume the Profile (Timer).
- **Stop key:** Press to stop any mode of operation.
- Page/Reset key: While in Setup Mode, press to access different Pages of configuration, Press this key to silence an alarm if the instrument alarm sounds during operation. In an operating mode, if an alarm or error condition occurs, press this key to return the instrument to normal operation once the condition is cleared.
- Menu/View key: While running any operating mode, pressing this key will display the high limit setpoint. While in Setup Mode, pressing this key will provide access to each Menu parameter.
- ▲▼ keys: Press these keys to adjust parameter settings. In Profile/Stopped Mode, press to select profile to run. In Profile/Run Mode, press ▲ key simultaneously with the Select key to force the program to advance one segment.

Outputs

The Protocol Plus control has seven different outputs available.

- Heating output: The control output is a DC voltage open-collector output which
 is time-proportioned and designed to control a heat control device such as a solid
 state relay.
- High limit: The high limit output is a form C relay which is energized under normal operating conditions. If the control senses a temperature higher than the high limit setpoint, or if there is a sensor error, the high limit relay will de-energize until the condition is cleared and the Reset key is pressed. When the high limit relay is de-energized, the heater is disabled.
- Retransmission: The retransmission output is a DC 1 to 5 volt or 4 to 20 ma
 (DC) signal that is proportional to the process temperature. The signal can be an
 input to other devices such as a chart recorder.
- Relay (four optional outputs): The four form A dry contact relay outputs can be configured to function as alarms, events, or end of cycle. These outputs can be utilized in Manual, Timer, or Profile Mode.

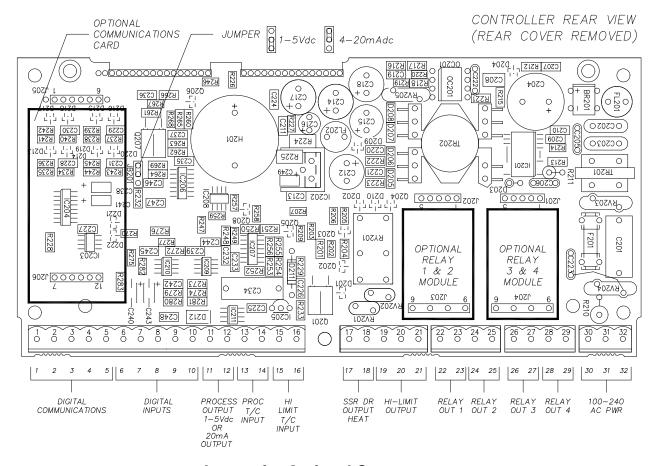

Layout for Optional Components

Relay (Continued)

Use the Relay Card Optional Ay p/n 144562 to add relays to the standard controller. Each relay card contains two relays (maximum of two cards Ay's allowed).

Communication

The Protocol Plus control has optional MODBUS communication available which can communicate via RS232, RS422, or RS485 to a computer. See communications option assembly p/n 161957 for board and cable assembly. Please refer to the MODBUS communications manual which comes with this option.

Optional Software

The Protocol Manager program allows the operator to start/stop multiple ovens (32 maximum) from a personal computer. A data log can also be used to record process information (p.n. 140008).

INSTRUCTIONS

Start-Up

These instructions are provided as a quick reference for operating the Protocol Plus control. If the Profile Mode is to be used, or the configuration of the control needs to be changed, please refer to the Setup Mode instructions before operating the control. For more detailed operating instructions refer to the Operation instructions for the mode you wish to use.

Upon initial power-up the control is in Manual/Stopped Mode (unless the Autostart or Fast Start Modes are active). To activate any operating mode from Stopped Mode, press the Select key until the desired mode is displayed, then press the Run key. If the proper Profile number is not displayed when the Profile Mode is accessed, press the ▲ or ▼ keys until the desired Profile number is displayed, then press the Run key. If no profile numbers can be displayed (display only reads NONE) then no profiles are currently programmed (see Setup Mode).

The Hi-limit thermocouple actual temperature reading is displayed, when the lower LCD display reads **HL Temp.** Note: This is not a error message.

The temperature setpoint can be adjusted while Manual or Timer Mode is running by pressing the UP or DOWN key.

To momentarily hold the Timer or Profile Mode, press the Hold key. To continue the Timer or Profile Mode, press the Run key.

To return to Stopped Mode at any time, press the Stop key and the cycle complete LED will illuminate.

Note that the control can be configured to automatically activate Manual, Timer or Profile Mode when power is applied (power switch turned on). See Control Page in the Setup Mode to utilize the Fast Start mode.

Operation

Manual Mode

Press the Select key until Manual is displayed (note you can press the Run key at any time to activate Manual Mode).

1. Press the Menu key to display the Process Temperature Setpoint (setpt). You can change the Setpoint with the ▲▼ keys.

Note: If the SPChange parameter on the Enable page in Setup Mode has been set to DISABLED, it must be changed to ENABLED before any changes to the process temperature and high limit setpoints can be made.

- Press the Menu key a second time to display current high limit setpoint (Hi-Lim SP). The value can be adjusted by pressing the ▲▼ keys. If Band is displayed, the high limit band feature is activated (see Control page) and the high limit can not be adjusted.
- 3. (optional feature) Press the Menu key a third time to display Event1. Press the ▲ key to turn on the event or ▼ to turn off the event. Repeat for all events which are enabled (up to 4).
- 4. To start Manual Mode, press the Run key.

The display will change from Stop to Run. To return to Stopped Mode, press the Stop key. While in operation, the process setpoint can be adjusted by using the keys to change the value while the mode is running. Pressing the Menu key will display the High Limit Setpoint (HLSP) setting.

If changes to the high limit setpoint or event output configuration are needed, they must be done from the stopped mode.

Timer Mode

- 1. Press the Select key until Timer is displayed (note you can press the Run key at any time to activate Timer Mode).
- Press the Menu key to display the Process Temperature Setpoint (Setpt). You can change the Setpoint with the ▲▼ keys.

Note that if the SPChange parameter on the Enable page in Setup Mode has been set to DISABLED, it must be changed to ENABLED before any changes to the process temperature and high limit setpoints can be made.

- 3. Press the Menu key a second time to display current high limit setpoint (Hi-lim SP). The value can be adjusted by pressing the ▲▼ keys. If **Band** is displayed, the high limit band feature is activated (see Control page) and the high limit can not be adjusted.
- 4. Press the Menu key a third time to display Time Set. You can change the time setting with the ▲▼ keys.
- 5. (optional feature) Press the Menu key a fourth time to display Event1. Press the ▲ key to turn on the event or ▼ to turn off the event. Repeat for all events which are enabled (up to 4).
- 6. Press the Menu key a fifth time to display the current guaranteed soak band (TmrGuarSoak) value. If the process temperature deviates from the setpoint by more than this value, the timer is placed in a hold condition. The timer continues when the process temperature falls within range. Reference the Quick Reference and Default Values section for available settings.
- 7. To start Timer Mode, press the Run key.

The display will change from Stop to Run and the time remaining will be displayed. To return to Stopped Mode, press the Stop key. While in operation, the process setpoint can be adjusted by using the $\blacktriangle \blacktriangledown$ keys to change the value while the mode is running. Pressing the Menu key will display the High Limit Setpoint.

Pressing the Run/Hold key while the Timer Mode is in operation will put the control in Hold status. The Timer LED will flash to indicate the held status. Press the Run/Hold key again to continue timing. The Timer LED will return to lit status.

Profile Mode

- 1. Press the Select key until Profile is displayed. "None" may be displayed if a profile has not been selected or no profiles entered.
- Press the ▲▼ key to display the desired profile to run.
- 3. To start Profile Mode, press the Run key.

The display will change from Stop to Run and the segment time remaining, Temperature Setpoint, Profile #, along with the current segment number, will be displayed. To return to Stopped Mode, press the Stop key.

Pressing the Run/Hold key while the Profile Mode is in operation will put the control in Hold status. Press the Run/Hold key again to continue the mode. The Profile LED will flash to indicate the hold status.

To advance to the next segment while running a profile, press the Select and UP arrow keys at the same time.

If **Link To** is set to **Standby** in the Program Page, at the End of Program/Profile,

- 1. Cycle Complete LED indication goes ON.
- 2. Controller beeps if **End of Cycle** beep is enabled.
- 3. Heater/control output keeps controlling oven temperature at last Soak setpoint.
- 4. All events programmed (if relay cards installed and programmed as an event) for the last Soak Segment stays active.

Note that ramping down too fast may cause the high limit relay to trip unexpectedly if the high limit band feature is used. This can be avoided by using a separate cooling profile that does not utilized the high limit band and then jumping to that profile to perform rapid cooling.

Auto Start Mode

The Auto Start Mode allows the control to start Manual, Timer, or Profile mode automatically at a preset time and day. See the Auto Start Page in Setup Mode for the time, day, and operating mode settings.

To activate the Auto Start Mode,

- 1. On Auto Start page, **Enable** is set to **Yes**.
- 2. LCD reads **Active** on line 1 in Auto Start Mode.
- 3. On Auto Start page **Enable** set to **No**, will deactivate Auto Start Mode.

Note that once you activate Auto Start, you can continue to use all operating modes as normal. If an operating mode is running at the time of a preset Auto Start function, and Auto Start is activated, the existing operating mode will override the auto Start function and the Auto Start will not turn on.

Note: All process Set to Run in Auto Start Mode must be at least one minute long for all Run Modes (Manual, Timer, and Profile).

Setup Mode

Configuration of the control and programming of the ramp/soak profiles must be done in the Setup Mode. To access Setup Mode, the control must first be in Stopped Mode.

- 1. Press the Select key until Setup is displayed.
- 2. Press the Page key and Security will be displayed.
- 3. Press the Menu key and Password will also be displayed. Use the ▲▼ keys to enter the proper password.
- 4. Once the proper password is displayed, press the Page key twice to enter the Setup Mode.

To exit Setup Mode, press and hold the Page key for three seconds.

The control has two levels of password-protected security. Level one provides access only to those menu pages that are enabled on the Enable page. Level two provides access to all menu pages, including the Enable page. The default security password values are 1 for level one and 2 for level two.

If an improper password has been entered, the control will remain at the Security display. To enter the proper password, press the Menu key. To exit Setup Mode, press and hold the Page key for three seconds.

Mapping of the Setup Mode is provided in the following sections. To access each parameter Page, which are described in detail in the following sections, press the Page key until the desired page heading is displayed. Press the Menu key to access each Menu parameter. Press the ▲▼ keys to change Menu parameter settings.

Refer to the Quick Reference and Default Values section for available settings for each Menu parameter.

Press the Page key to continue with each Page, or press and hold the Page key for three seconds to exit Setup Mode.

Instructions for Setup Mode Pages

Program Page

Programming of the profiles is provided on the Program Page. Eight profiles are available with up to eight ramp and soak segments per profile.

If the optional relay outputs are installed, they must be configured as alarms or events on the Relay Outputs Page before they can be utilized. If configured as event outputs, these relays can be used as time or temperature events.

When entering the Program Page, press the Select key to select the profile you wish to enter/edit, then press the Menu key. The first parameter (Profile #, Segment 1, Ramp Time) will display. Adjust the time value with the ▲▼ keys. Once the proper value is displayed, press the Menu key to continue. Continue with the Menu key to adjust/view each parameter.

If the ramp time value of the current segment is left at 0:00, the next press of the Menu key will advance the control to the High Limit Setpoint parameter for that profile. Continue entering / verifying all parameters until you get to the last parameter (Guaranteed Soak Band). Once all parameters have been properly entered, press the Page key to return to the top of the Profile Page. You can press the Select key to enter/edit another profile, press the Page key to access another page, or press and hold the Page key to exit Setup mode.

While editing any profile, pressing the Select key will advance the control to the time value for the next segment, until the last segment has been reached. This allows faster editing of the profile rather than pressing the Menu key to advance past each parameter.

If **Link To** is set to **Standby** in the Program Page, at the End of Program/Profile,

- 1. Cycle Complete LED indication goes ON.
- 2. Controller beeps if **End of Cycle** beep is enabled.
- 3. Heater/control output keeps controlling oven temperature at last Soak setpoint.
- 4. All events programmed (if relay cards installed and programmed as an event) for the last Soak Segment stays active.

To run a profile indefinitely, link the profile to itself.

Menu Item	Display	Description
Ramp Time Seg 1	Pro-1 Seg-1 Ramp Time	Ramp time for segment 1 of profile
Event 1 Set Value*	Pro-1 Seg-1 Ramp Event 1	Event 1 setting for segment 1 ramp of profile
Event 2 Set Value*	Pro-1 Seg-1 Ramp Event 2	Event 2 setting for segment 1 ramp of profile
Event 3 Set Value*	Pro-1 Seg-1 Ramp Event 3	Event 3 setting for segment 1 ramp of profile
Event 4 Set Value*	Pro-1 Seg-1 Ramp Event 4	Event 4 setting for segment 1 ramp of profile
Soak Temp Seg 1	Pro-1 Seg 1 Soak Temp	Soak temperature for segment 1 of profile
Soak Time Seg 1	Pro-1 Seg 1 Soak Time	Soak time for segment 1 of profile
Event 1 Set Value*	Pro-1 Seg-1 Soak Event 1	Event 1 setting for segment 1 soak of profile
Event 2 Set Value*	Pro-1 Seg-1 Soak Event 2	Event 2 setting for segment 1 soak of profile
Event 3 Set Value*	Pro-1 Seg-1 Soak Event 3	Event 3 setting for segment 1 soak of profile
Event 4 Set Value*	Pro-1 Seg-1 Soak Event 4	Event 4 setting for segment 1 soak of profile
(repeat for segments 2	2-8, until ramp or soak time = 00:00	0)
High Limit Setpoint	Pro-1 Hi-Lim SP	High limit setpoint for profile**
Loop From	Pro-1 Loop From Seg	To start a loop action in a profile
Loop To	Pro-1 Loop To Seg	To end a loop action in a profile
Loop Count	Pro-1 Loop Number	Number of times to execute loop
Profile Link	Pro-1 Link To Pro	To jump from this profile to another

See the definitions on the following pages for parameter ranges.

Guaranteed soak band for profile

Guaranteed Soak

Pro-1 Guar Band

^{*} only available if optional relay outputs are installed (repeat all for profiles 2-8)

^{**} Set to **Band** to use the high limit band feature

There are eight profiles available. Profile #

Segment# Recipe segments 1 through 8 may be programmed, each with its own

set of events, ramp and soak times, and soak temperature.

Ramp Time The time required to ramp from one setpoint up to another setpoint.

> Values between 0 and 99:59 are allowable. In the Protocol Plus controller, the profile ramp and soak times are stored without units. Units are set as either hours and minutes (HH:MM) or minutes and seconds (MM:SS). The setpoint will automatically increment from the

actual temperature to the soak temperature.

EV1 through 4 From 1 to 4 events may be programmed into the ramp time portion of

> each segment here. These typically involve actuating/disabling relays to close/open valves or perform other relay-controlled functions.

> NOTE: These will only actuate when the controller has the relay cards

installed and programmed for an event.

Soak Temp. The temperature setpoint of a particular segment is entered here; it

can range from -18 to 540 degrees C (0 to 1000 degrees F).

Soak Time The duration of soak is entered here; the value can range from 0 to

99:59.

EV1 through 4 From 1 to 4 events may be programmed into the soak portion of each

> segment here. These typically involve actuating/disabling relays to close/open valves or perform other relay-controlled functions. NOTE:

These will only actuate when the controller has the relay cards

installed and programmed for an event.

Hi Limit SP The high limit setpoint may be entered here; if the temperature

exceeds this value, the hi-limit will alarm and shut off the heater.

Loop From Values are No, Seg-1 to Seg-8.

Loop To Values are No, Seg-1 to Seg-8.

Loop Number Values are 0 - 99.

These values enable the operator to jump from a certain step to

another step of the recipe a preset number of times.

Profile Link Values are STANDBY/STOP/HOLD/1 - 8. When the profile ends, the

> profile can hold the temperature setpoint while keeping the events active, turn the heater off, hold the temperature setpoint at the end of


the profile, or jump to another specified profile.

Guaranteed If the process temperature deviates from the setpoint by more than Soak Band

this value, the soak timer is placed in a hold condition. The timer

continues when the process temperature falls within range.

Sample Profile

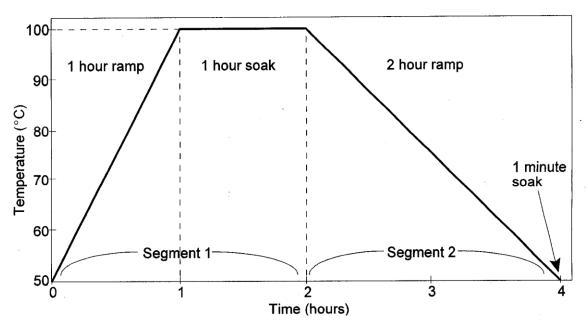

Programming Table

Profile	Number	_1		Pro	ofile Na	me					
Seg-		F	Ramp					Soak			
ment	Time	1		ents		Temp-	Time			ents	
		1	2	3	4	erature		1	2	3	4
1	01h00					100	01h00				
2	02h00					50	00h01				
3	00h00										
4											
5											
6											
7											
8											
	High Limit	Setpoint		115							
	Loop Froi	m Seg		No							
	Loop To	Seg		No							
	Loop Nu	mber		0							
	Link To	Го Рго									
Guar Soak Band 10											

Auto Start Page

The Auto Start Page can be configured to automatically start Manual, Timer or Profile Mode at a specified time and day. Note that if Auto Start Enable is set to Yes in the Setup Mode, the Auto Start feature is not turned on - it is available to the operator to be activated in Stopped Mode.

To configure the Auto Start feature:

- 1. Access the Setup Mode.
- Press the Page key until Auto Start is displayed.
- 3. Press the Menu key. If there is no change in the display, the controller may not have the realtime clock option.
- 4. Set Auto Start Enable to Yes.
- 5. Using the Menu key, scroll through the options available and use the ▲▼ keys to set the mode desired for each day of the week. You may select from Manual, Timer, or Profile 1 through 8.
- 6. When the mode is set press the Menu key.
- 7. Enter the time of day you wish the mode to activate.
- 8. Continue through the rest of the week by pressing the Menu key, or press the Page key when done.

One Auto Start mode can be set for each day of the week. Exit the Setup mode by pressing and holding the Page key for three seconds. Press the Select key until Auto Start is displayed. Make sure the correct time and day is displayed. If not proper, set it to the correct time on the Real Time Clock Page in the Setup mode.

To activate the Auto Start Mode,

- 1. On Auto Start page, **Enable** is set to **Yes**.
- 2. LCD reads Active on Line 1 in Auto Start Mode.
- 3. On Auto Start page **Enable** set to **No**, will deactivate Auto Start Mode.

Note that once you activate Auto Start, you can continue to use all operating modes as normal. If an operating mode is running at the time of a preset Auto Start function, and Auto Start is activated, the existing operating mode will override the Auto Start function and the Auto Start will not turn on.

Note: All process Set to Run in Auto Start Mode must be at least one minute long for all Run Modes (Manual, Timer, and Profile).

Menu Item	Display	Description	Range
Enable Autostart	Auto Start Enable	Enable (yes) or disable (no) the Autostart function	No, Yes
Sunday mode	Auto Start Sun Mode	Set mode on Sunday to activate	Off, Manual, Timer, Pro-1 to Pro-8
Sunday time	Auto Start Sun Time	Set time on Sunday for mode to activate	00:00 to 23:59
Monday mode	Auto Start Mon Mode	Set mode on Monday to activate	Off, Manual, Timer, Pro-1 to Pro-8
Monday time	Auto Start Mon Time	Set time on Monday for mode to activate	00:00 to 23:59
Tuesday mode	Auto Start Tue Mode	Set mode on Tuesday to activate	Off, Manual, Timer, Pro-1 to Pro-8
Tuesday time	Auto Start Tue Time	Set time on Tuesday for mode to activate	00:00 to 23:59
Wednesday mode	Auto Start Wed Mode	Set mode on Wednesday to activate	Off, Manual, Timer, Pro-1 to Pro-8
Wednesday time	Auto Start Wed Time	Set time on Wednesday for mode to activate	00:00 to 23:59
Thursday mode	Auto Start Thu Mode	Set mode on Thursday to activate	Off, Manual, Timer, Pro-1 to Pro-8
Thursday time	Auto Start Thu Time	Set time on Thursday for mode to activate	00:00 to 23:59
Friday mode	Auto Start Fri Mode	Set mode on Friday to activate	Off, Manual, Timer, Pro-1 to Pro-8
Friday time	Auto Start Fri Time	Set time on Friday for mode to activate	00:00 to 23:59
Saturday mode	Auto Start Sat Mode	Set mode on Saturday to activate	Off, Manual, Timer, Pro-1 to Pro-8
Saturday time	Auto Start Sat Time	Set time on Saturday for mode to activate	00:00 to 23:59

PID Page

The PID Page contains parameters which control the response to the setpoint and process variable input. To access the PID Page, enter the Setup Mode. Press the Page key until PID is displayed. Press the Menu key. Each parameter can be changed by pressing the Menu key until the desired parameter is displayed, and then pressing the $\blacktriangle \nabla$ keys to change the value.

Menu Item	Display	Description	Range
Display units	PID Temp Unit	Set display units to °C or °F	°C or °F
Proportional band	PID Prop Band	Set proportional band for tuning	1 to 56°C (1 to 100°F)
Integral reset	PID Reset/Rep/Min	Set integral reset for tuning	0.0 to 100 seconds/repeat
Derivative Rate	PID Rate In Sec	Set derivative rate for tuning	0.0 to 500 seconds
AutoTune	PID AutoTune	Enable auto tuning function	Disable, Enable

The AutoTune parameter disables or enables the AutoTune function. To utilize AutoTuning:

- 1. Access the Setup Mode.
- Press the Page key until the display reads AutoTune. Press the ▲ key to enable the AutoTune.
- 3. Press the Page key for three seconds to exit Setup Mode.
- 4. Cycle power to the instrument.
- 5. Set Manual Mode to run. The display will alternately display AutoTune and Manual.

If the Manual Mode setpoint is less than 50 degrees higher than the actual process temperature, the AutoTune function will create an error condition. This can be cleared by either cooling off the process temperature or increasing the setpoint until there is more than 50 degrees between them. Once the AutoTune function is allowed to complete tuning, the AutoTune parameter will disable by itself.

If you wish to cancel the AutoTune function, press the STOP key, access the PID page in Setup Mode, and set the AutoTune parameter to Disable.

Control Page

The Control Page contains various parameters which control miscellaneous functions. To access the Control Page, enter the Setup Mode. Press the Page key until Control is displayed. Press the Menu key. Each parameter can be changed by pressing the Menu key until the desired parameter is displayed, and then pressing the ▲▼ keys to change the value.

Menu Item	Display	Description	Range
Cycle Time	Control Cycle Time Sec	Set cycle time in seconds for control output	1 to 60 seconds
High limit setpoint	Control Hi-Lim SP***	Maximum value for all high limit setpoints	MinHiLimSP - MaxHiLimSP*
High limit band	Control Hi-Lim Band	If=0, high limit setpoint= Control Hi-Lim SP If>0, high limit setpoint= Control SP* + Band	Off, 3°C to 11°C (5°F to 20°F)
Power fail recovery	Control PwrFRec	Controls response to loss of power	Stop, Restart, Hold, Resume
Recovery time limit	Control PFRTime	Control aborts to Stopped mode if power is lost for time period longer then set value	00m00s to 99m59s
Powerup start enable	ControlPwrUpStrt	Allows mode to automatically start when power is first applied	Disable, Enable
Powerup Start Mode	Control StrtMode	Operating mode for powerup start	Off, Manual, Timer, Pro-1 to Pro-8
Hysteresis	Control Hysteresis	Hysteresis for all alarms and temperature events	1°C to 56°C (1°F to 100°F)
Process out low	Control RetOutLo	Process value for retransmit output = 1VDC	-73°C to 760°C (-100°F to 1400°F)
Process out high	Control RetOutHi	Process value for retransmit output = 5VDC	-73°C to 760°C (-100°F to 1400°F)
Time scale	Control TimeScale	Time scale setting for program and timer mode**	hh:mm or mm:ss
Key press beep	Control KeyBeep	Internal beeper sounds when key is pressed	On or Off
End of cycle beep	Control EOCBeep	Internal beeper sounds at end of cycle	On or Off
Alarm beep	Control AlarmBeep	Internal beeper sounds for alarms	On or Off

^{*} includes ramping setpoints during profiles and controlled ramps

^{**} power fail recovery time limit is always MM:SS regardless of time scale setting

^{***} high limit setpoint is a read-only item which is calculated on Enable page

Communication Page (optional)

The Communication Page contains parameters for the communications feature. To access the Communications Page, enter the Setup Mode (see description earlier in this manual). Press the Page key until Communication is displayed. Press the Menu key. (NOTE: If there is no change in the display, the controller does not have the communications board installed.) Each parameter can be changed by pressing the Menu key until the desired parameter is displayed, and then pressing the $\blacktriangle \nabla$ keys to change the value.

Menu Item	Display	Description	Range
Address	Communication CommAddr	Sets address node for control	1 to 247
Mode	Communication Mode	Turns on/off communications	OFF, Modbus
Baud Rate	Communication BaudRate	Sets interface speed	2400, 4800, 9600, 19.2K, 38.4K
Parity	Communication Parity	Sets parity for interface	None, Odd, Even

Real Time Clock Page

The Real Time Clock Page allows the control to be configured to have an operating mode begin automatically at a specific time on a specific day of the week. The real time clock feature is also used for using the Power Failure Recovery mode Time Limit feature (see Control Page). The real time clock is a seven day, 24 hour clock with battery backup.

To access the Real Time Clock Page, enter the Setup Mode. Press the Page key until Clock is displayed. Press the Menu key. Each parameter can be changed by pressing the Menu key until the desired parameter is displayed, and then pressing the ▲▼ keys to change the value.

Menu Item	Display	Description	Range
Day of the week	Clock Day	Setting for current day of the week	Sun, Mon, Tue, Wed, Thu, Fri, Sat
Time of day	Clock HH:MM	Setting for current time of the day	00:00 to 23:59
Reset clock	Clock UP to Reset CLK*	Press the key to set the clock to entered values	Ready, Done

^{*} If the \triangle key is not pressed, the clock values will retain their original values. The display will change to Done if the clock is reset

Relay Outputs Page (optional)

The Relay Outputs Page configures the four alarm/event outputs. Each output has a dedicated indicator light and can be configured as a temperature alarm, profile event, or end of cycle output. Temperature alarms can be of type process high, process low, deviation high, deviation low, or deviation band.

To access the Relay Page, enter the Setup Mode (see description earlier in this manual). Press the Page key until Relay is displayed. Press the Select key until the desired relay output is selected. Press the Menu key. Each parameter can be changed by pressing the Menu key until the desired parameter is displayed, and then pressing the ▲▼ keys to change the value. To configure a specific relay, press the Select key until the desired relay appears.

NOTE: If Relay 0 appears, then no relays are installed (see relay kit assembly p.n. 144562).

Menu Item	Display	Description	Range
Type of relay	Relay 1	Set function of relay	Off, Alarm, Cycl, Ev1 to Ev4
	RelayType		
Action of relay	Relay 1	Set coil and contact state of	NDE, NE, NDEL, NEL****
	RelayAction	relay	
Type of	Relay 1	Set alarm type for relay	High, Low, Plus, Minus, Band
alarm*	AlarmType	, , , , , , , , , , , , , , , , , , ,	
Alarm	Relay 1 AlmHi/Lo	Setpoint for alarm	-73°C to 760°C (-100°F to 1400°F)
setpoint*	SP	·	
Alarm	Relay 1	Deviation band for alarm	1 to 56°C (1 to 100°F)
deviation*	AlmDevBand		,
Inhibit alarm*	Relay 1	Inhibits alarm until "safe"	En or Dis
	ALMInhibit	condition is reached	
Type of	Relay 1	Set event type for relay	Time or Temp
event**	EventType		
Event	Relay 1 Event SP	Setpoint for temperature event	SPLoLim to SPUpLim****
setpoint***			·

(repeat for relay outputs 2-4, if available)

Turning on the Alarm Inhibit function disables the alarm output on power up until the process temperature has reached a non-alarming condition ("safe").

If the relay output has been configured as latching, the RESET key must be pressed to return the output to the non-alarm state once the alarm condition has cleared.

^{*} appears only for alarm types

^{**} appears only for time or temperature event types

^{***} appears only for temperature event types

Test Page

The Test Page contains parameters which allow manual control of the heat control and optional relay outputs and should be used only for testing the functionality of the control instrument. Do not operate the oven for processes using the Test Page.

To access the Test Page, enter the Setup Mode (see description earlier in this manual). Press the Page key until Test is displayed. Press the Menu key. Each parameter can be changed by pressing the Menu key until the desired parameter is displayed, and then pressing the $\blacktriangle \blacktriangledown$ keys to change the value.

Menu Item	Display	Description	Range
Heater	Test HeatOut	Activate SSR output 100%	On
output			
High limit	Test HiLimOut	Activate high limit alarm	On
relay		(de-energize relay)	
Relay 1	Test Rly1 Out	Energize relay output 1	On
output			
Relay 2	Test Rly2 Out	Energize relay output 2	On
output			
Relay 3	Test Rly3 Out	Energize relay output 3	On
output			
Relay 4	Test Rly4 Out	Energize relay output 4	On
output			
HiLim	Test HL Temp	Displays sensor reading*	
Sensor	(push and hold up)		

^{*}Push ▲ key to refresh display

When the Test Page is entered, all outputs are automatically set to off. When exiting the Test Page, all outputs will return to their previous condition regardless of the Test Page settings.

Zone Calibration Page

The Zone Calibration Page allows adjustment of the displayed temperature versus the actual temperature measured by the control thermocouple. This may be desirable in certain conditions where the center of the oven chamber is not the same temperature as the control thermocouple. This may occur when the oven is not allowed to soak at a constant temperature for long periods of time, or the oven is being used at high temperature.

There is also a Factory Calibration Recovery which will restore the factory calibration values when the control was first shipped by the manufacturer. This may be helpful if the calibration has been lost and a calibration instrument is not readily available. To use the Factory Calibration Recovery feature only, bypass the Zone 1 and Zone 2 calibration parameters by pressing the Menu key.

To access the Zone Calibration Page, enter the Setup Mode (see description earlier in this manual). Press the Page key until Zone Cal is displayed. Press the Menu key. Each parameter can be changed by pressing the Menu key until the desired parameter is displayed, and then pressing the $\blacktriangle \blacktriangledown$ keys to change the value.

Menu Item	Display	Description	Range
Zone 1	Zone Cal	Point at which Zone 1 is set (center	-73°C to 760°C (-100°F to 1400°F)
actual	Zone1Act	of chamber)	
Zone1	Zone Cal	Desired displayed value for Zone 1	-73°C to 760°C (-100°F to 1400°F)
displayed	Zone1Dis	setting	
Zone 2	Zone Cal	Point at which Zone 2 is set (center	-73°C to 760°C (-100°F to 1400°F)
actual	Zone2Act	of chamber)	
Zone2	Zone Cal	Desired displayed value for Zone 2	-73°C to 760°C (-100°F to 1400°F)
displayed	Zone2Dis	setting	
Factory	Zone Cal	Restores the factory calibration	Ready or Done (push- key)
calibration*	FactCal	values	

^{*}Only use when no calibration instrument is available, push ▲ key to restore factory values.

Press the Page key to exit the Zone Calibration Page.

Two points of display calibration (temperature offset) are available. The Zone 1 Actual and Zone 2 Actual parameters are the two temperature points where the offset is to take effect. These values are adjustable. The Zone 1 and Zone 2 Displayed parameters are the values the user wishes to have displayed at the Actual temperatures, and are also adjustable.

As an example, the control is displaying 400°F with the setpoint being 400°F, but the center of the oven chamber is actually 395°F. This can occur due to oven wall losses and product loading variations. The operator may change the zone calibration so that the center of the oven is 400°F when the display reads 400°F. In this case operate the oven in manual mode with a setpoint of 400°F. Record the center of the chamber (as measured with an independent sensor). Access the Zone Calibration Page and enter this measured value as the Zone 2 Actual value, with 400° as the Zone 2 Displayed value.

Zone 1 can be adjusted using the same method at a lower temperature. The instrument will then create a linear offset based on the Zone1 and Zone 2 Actual temperature values. Note that the oven does not have to be heated to adjust the zone parameters if the zone values are known based on prior experience.

Sensor Calibration Page

The Sensor Calibration Page has parameters which can change the internal calibration of the temperature sensor input signal. There is a low and high calibration point for both the control sensor and the high limit sensor. To calibrate the instrument, allow the control to warm up for at least 30 minutes.

To access the Sensor Calibration Page, enter the Setup Mode (see description earlier in this manual). Press the Page key until Control Sensor is displayed. Press the Menu key.

The control may have the optional process value retransmission output feature. The output is a 1 to 5VDC signal. To calibrate the retransmit signal, the RetOutLo and RetOutHi values from the Control Page must be known. You may bypass calibrating the control and high limit sensor input to access only the retransmit calibration by pressing the Menu key until RetCalLo is displayed (skip steps 4-17).

To re-calibrate the instrument:

- 1. Disconnect the control and high limit sensor thermocouples.
- 2. Connect a calibration instrument with a type J thermocouple output to the control sensor input. Allow the control to warm up at least 30 minutes.
- 3. Access Setup Mode.
- 4. Press Page key until -100F is displayed.
- 5. Press Menu key until **Ctrl Sens -100F** is displayed.
- 6. With **Ctrl Sens -100F** displayed, adjust the calibration instrument to Type J thermocouple, -100 degrees Fahrenheit output.
- 7. Wait 30 seconds. Press the ▲ key.
- 8. With **Ctrl Sens 1400F** displayed, adjust the calibration instrument to 1400 degrees Fahrenheit output (Type J thermocouple).
- 9. Wait 30 seconds. Press the ▲ key.
- 10. When the control displays **Ctrl Sens Done**, disconnect the calibration instrument from the control sensor input and connect it to the high limit sensor input. Reconnect the control sensor thermocouple.
- 11. Press the Menu key until **HL Sens** -100**F** is displayed.
- 12. With the control displaying **HL Sens -100F**, adjust the calibration instrument to -100 degrees Fahrenheit output (Type J thermocouple).
- 13. Wait 30 seconds. Press the ▲ key.
- 14. With the control displaying **HL Sens 1400F**, adjust the calibration instrument to 1400 degrees Fahrenheit output (Type J thermocouple).
- 15. Wait 30 seconds. Press the ▲ key.
- When the control displays HL Sens Done, disconnect the calibration instrument from the high limit sensor input. Re-connect the high limit sensor thermocouple.
- 17. To skip calibration of the retransmit signal, press the Page key twice to exit the Sensor Calibration Page.
- 18. To calibrate the retransmit signal, press the Menu key until **RetCalLo** is displayed.

- 19. Connect a calibration instrument with a type J thermocouple output to the control sensor input.
- 20. Connect a voltage measurement device to the retransmit output terminals.
- 21. Set the calibration instrument output to the temperature value of the **RetOutLo** parameter from the Control Page.
- 22. Adjust the **RetCalLo** * value using the ▲▼ keys until the voltage measurement device reads 1VDC.
- 23. Press the Menu key.
- 24. Set the calibration instrument output to the temperature value of the **RetOutHi** parameter from the Control Page.
- 25. Adjust the **RetCalHi** * value using the ▲▼ keys until the voltage measurement device reads 5VDC.
- 26. Press the Menu key.
- 27. Press the Page key to exit the Sensor Calibration Page.
- 28. Calibration is now complete. Disconnect the calibration instrument and voltage measurement device (if used).
- 29. Verify that the control and high limit sensor thermocouples are connected.

Menu Item	Display	Description	Range
Control Sensor	Ctrl Sens -100F	Calibrate Sensor Low End	-100 to 1400°F
Cal	Ctrl Sens 1400F	Calibrate Sensor High End	-100 to 1400°F
	Ctrl Sens Done	Control Sensor Cal Complete	(read only)
HiLim Sensor	HL Sens -100F	Calibrate HiLim Sensor Low	-100 to 1400°F
Cal	HL Sens 1400F	End	-100 to 1400°F
	HL Sens Done	Calibrate HiLim Sensor High	(read only)
		End	
		HiLim Sensor Cal Complete	
Retransmit Cal	RetCalLo XXXX *	Calibrate Retransmit Output	0 to 4096**
	RetCalHi XXXX *	Low	0 to 4096***
		Calibrate Retransmit Output	
		High	

^{*}Note that the actual RetCalLo and RetCalHi values displayed are of no importance.

Enable Page

The Enable Page controls access to the other Setup Pages. The setpoint minimum and maximum values, and the security passwords are also set on the Enable Page. To access the Enable Page, enter the Setup Mode using a level 2 access code (see description earlier in this manual). Press the Page key until Enable is displayed. Press the Menu key. Each parameter can be changed by pressing the Menu key until the desired parameter is displayed, and then pressing the AV keys to change the value.

NOTE: Changing the enable to "yes" for each page will allow access to the page in Level 1 security.

Menu Item	Display	Description	Range
Profiles	Enable Profile 1-8	Controls access to Program Page	Yes or No
Autostart	Enable Auto Start	Controls access to AutoStart Page	Yes or No **
PID	Enable PID	Controls access to PID Page	Yes or No
Control	Enable Control	Controls access to Control Page	Yes or No
Communication	Enable Communication	Controls access to Communication Page	Yes or No **
Real Time Clock	Enable Clock	Controls access to Real Time Clock Page	Yes or No **
Relay outputs	Enable Relay 1-4	Controls access to Relay Page	Yes or No **
Test	Enable Test	Controls access to Test Page	Yes or No
Zone Calibration	Enable Zone Cal	Controls access to Zone Calibration Page	Yes or No
Sensor Calibration	Enable Sensor Cal	Controls access to Sensor Calibration Page	Yes or No
Setpoint lower limit	Enable SPLowerLim	Sets minimum setpoint allowed	-73°C to 759°C (-100°F to 1399°F)
Setpoint upper limit	Enable SPUpperLim	Sets maximum setpoint allowed	-73°C to 760°C (-100°F to 1400°F)
High limit overhead	Enable HiLimOH	Sets maximum high limit setpoint allowed*	3 to 11°C (5 to 20°F)*
Password level	Enable Password 1	Sets password for access level 1	0 to 999
Password level 2	Enable Password 2	Sets password for access level 2	0 to 999
Setpoint Change	Enable SPChange	Set to DISABLE to lock out setpoint and high limit setpoint changes in Manual and Timer Modes	Yes or No
Analog Output Type	Enable Analog Type	Sets Analog Output type	Ctrl or Proc

^{*} Maximum high limit setpoint = SPUpperLim + HiLimOH

^{**} If the controller does not have this feature, "No" is the only option.

Digital Inputs (optional)


The Protocol Plus control can be run by external inputs wired to the control from an external source such as a PLC or control panel switches. The external run operation can Run, Hold or Stop profiles 1 through 7 (profile 8 can not be operated externally). Refer to the table below for the inputs required for the desired operation. NOTE: A profile must be created in the program page before trying to run a profile number.

Input 1	Input 2	Input 3	Profile Selected
ON	OFF	OFF	1
OFF	ON	OFF	2
ON	ON	OFF	3
OFF	OFF	ON	4
ON	OFF	ON	5
OFF	ON	ON	6
ON	ON	ON	7
OFF	OFF	OFF	none

To start the selected profile, set Input 1, 2, 3 or 4 to ON.

To hold a profile, set Input 4 to OFF.

To stop a profile, set all inputs to OFF.

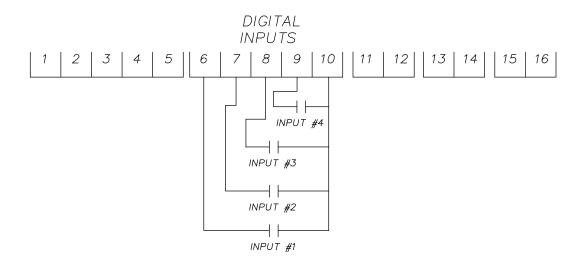

Digital Inputs

Error Messages and Alarms

The Alarm Status <u>Hi-limit</u> LED is flashing. This indicates a problem with the thermocouple, or the Hi-limit setpoint has been exceeded. Once the problem has corrected, press the Reset pushbutton.

The Alarm Status **Soak** LED is flashing. This indicates that the oven temperature has not entered or dropped out of the soak band and the soak timer has stopped.

The top LED Display reads **OPEN** and the lower LCD display reads **CONTROL SENS ERR**. This indicates that the Control thermocouple is disconnected or broken. Repair or replace the thermocouple.

The lower LCD display reads **HI-LIM SENS ERR**. This indicates that the Hi-limit thermocouple is disconnected or broken. Repair or replace the thermocouple.

The lower LCD display reads **HIGH LIMIT ALARM**. This indicates that the Hi-limit temperature setpoint has been exceeded. Determine if the setting is set too close to the setpoint, the SSR is defective, or the calibration is incorrect.

Note: The lower LCD display intermittently reads **HL Temp.** This is not an error message, but the Hi-limit thermocouple temperature reading.

Quick Reference and Default Values

Program Page

Menu Item	Display	Default	Range	Setting
Ramp Time Seg 1	Pro-1 Seg-1 Ramp Time	00:00	00m00s to 99h59s	
Event 1 Set Value	Pro-1 Seg-1 Ramp Event 1	Off	Off, On	
Event 2 Set Value	Pro-1 Seg-1 Ramp Event 2	Off	Off, On	
Event 3 Set Value	Pro-1 Seg-1 Ramp Event 3	Off	Off, On	
Event 4 Set Value	Pro-1 Seg-1 Ramp Event 4	Off	Off, On	
Soak Temp Seg 1	Pro-1 Seg 1 Soak Temp	68°F	SPLowerLim to SPUpperLim *	
Soak Time Seg 1	Pro-1 Seg 1 Soak Time	00:00	00m00s to 99h59s	
Event 1 Set Value	Pro-1 Seg-1 Soak Event 1	Off	Off, On	
Event 2 Set Value	Pro-1 Seg-1 Soak Event 2	Off	Off, On	
Event 3 Set Value	Pro-1 Seg-1 Soak Event 3	Off	Off, On	
Event 4 Set Value	Pro-1 Seg-1 Soak Event 4	Off	Off, On	
(repeat for segments	2-8)		· I	
High Limit Setpoint	Pro-1 Hi-Lim SP	Max HiLimSP	MinHiLimSP to MaxHiLimSP *, Band **	
Loop From	Pro-1 Loop From XX	No	No, Seg-1 to Seg-8	
Loop To	Pro-1 Loop To XX	No	No, Seg-1 to Seg-8	
Loop Count	Pro-1 Loop Number	0	0 to 99	
Profile Link	Pro-1 Link To XX	Stop	Standby***, Stop, Hold, Pro-1 to Pro-8	
Guaranteed Soak	Pro-1 Guar Band	Off	Off, 1 to 778°C	

^{*} See Enable Page

- 1. Cycle Complete LED indication goes ON.
- 2. Controller beeps if End Of Cycle beep is enabled.
- 3. Heater/control output keeps controlling oven temperature at last Soak setpoint.
- 4. All events programmed (if relay cards installed and programmed as an event) for the last Soak Segment stays active.

^{**} Band value is set on Control Page

^{***} If Standby

Programming Table

Profile	Number	

Profile Name_____

Seg-		F	Ramp					Soak			
ment	Time		Eve	ents		Temp-	Time			ents	
		1	2	3	4	erature		1	2	3	4
1											
2											
3											
4											
5											
6											
7											
8											
	High Limit S	Setpoint									
	Loop Fror	m Seg									
	Loop To	Seg									
	Loop Nu	mber									
	Link To	Pro									
	Guar Soal	k Band									

Autostart

Menu Item	Display	Default	Range	Setting
Enable Autostart	Auto Start Enable	No	No, Yes	
Sunday mode	Auto Start Sun Mode	Off	Off, Manual, Timer, Pro-1 to Pro-8	
Sunday time	Auto Start Sun Time	00:00	00:00 to 23:59	
Monday mode	Auto Start Mon Mode	Off	Off, Manual, Timer, Pro-1 to Pro-8	
Monday time	Auto Start Mon Time	00:00	00:00 to 23:59	
Tuesday mode	Auto Start Tue Mode	Off	Off, Manual, Timer, Pro-1 to Pro-8	
Tuesday time	Auto Start Tue Time	00:00	00:00 to 23:59	
Wednesday mode	Auto Start Wed Mode	Off	Off, Manual, Timer, Pro-1 to Pro-8	
Wednesday time	Auto Start Wed Time	00:00	00:00 to 23:59	
Thursday mode	Auto Start Thu Mode	Off	Off, Manual, Timer, Pro-1 to Pro-8	
Thursday time	Auto Start Thu Time	00:00	00:00 to 23:59	
Friday mode	Auto Start Fri Mode	Off	Off, Manual, Timer, Pro-1 to Pro-8	
Friday time	Auto Start Fri Time	00:00	00:00 to 23:59	
Saturday mode	Auto Start Sat Mode	Off	Off, Manual, Timer, Pro-1 to Pro-8	
Saturday time	Auto Start Sat Time	00:00	00:00 to 23:59	

PID

: ·-				·
Menu Item	Display	Default	Range	Setting
Display units	PID Temp Unit	°C	°C or °F	
Proportional band	PID Prop Band	6°C	1 to 56°C (1 to 100°F)	
Integral reset	PID Reset	2	0 to 100 seconds/repeat	
Derivative Rate	PID Rate	0	0 to 500 seconds	
AutoTune	PID AutoTune	Disable	Disable, Enable	

Control

Menu Item	Display	Default	Range	Setting
Cycle Time	Control Cycle Time	1	1 to 60 seconds	
High limit setpoint	Control Hi-Lim SP	Max HiLimSP	MinHiLimSP - MaxHiLimSP*	
High limit band	Control Hi-Lim Band	Off	Off, 3°C to 11°C (5°F to 20°F)	
Power fail recovery	Control PwrFRec	Stop	Stop, Restart, Hold, Resume	
Recovery time limit	Control PwrFTime	00m00s	00m00s to 99m59s	
Powerup start enable	Control EPwrStrt	Dis	Dis, En	
Powerup Start Mode	Control StrtMode	Off	Off, Manual, Timer, Pro-1 to Pro-8	
Hysteresis	Control Hyst	3°C	1°C to 56°C (1°F to 100°F)	
Process out low	Control RetOutLo	80°C	-73°C to 760°C (-100°F to 1400°F)	
Process out high	Control RetOutHi	400°C	-73°C to 760°C (-100°F to 1400°F)	
Time scale	Control TimeScale	hh:mm	hh:mm or mm:ss	
Key press beep	Control KeyBeep	On	On or Off	
End of cycle beep	Control EOCBeep	Off	On or Off	
Alarm beep	Control AlarmBeep	Off	On or Off	

*see Enable Page

Communication (optional)

Communication	optional)			
Menu Item	Display	Default	Range	Setting
Address	Communication CommAddr	1	1 to 247	
Mode	Communication CommMode	OFF	OFF, Modbus	
Baud rate	Communication Baud Rate	19.2K	2400, 4800, 9600, 19.2K, 38.4K	
Parity	Communication Parity	None	None, Odd, Even	

Real Time Clock

Menu Item	Display	Default	Range	Setting
Day of the week	Clock Day	Mon	Sun, Mon, Tue,	-
			Wed, Thu, Fri, Sat	
Time of day	Clock HH:MM	00:00	00:00 to 23:59	
·				
Reset clock	Clock UP to Reset CLK*	Ready	Ready, Done	

^{*} if the \triangle key is not pressed, the clock values will retain their original values, the display will change to Done if the clock is reset

Relay Outputs (optional)

Push Select key to select relay. If Relay 0 appears, no relays are installed

Menu Item	Display	Default	Range
Type of relay	Relay 1 RelayType	Off	Off, Alarm, Cycl, Ev1 to Ev4
Action of relay	Relay 1 RelayAction	NDE	NDE, NE, NDEL, NEL****
Type of alarm*	Relay 1 AlarmType	High	High, Low, Plus, Minus, Band
Alarm setpoint*	Relay 1 AlmHi/Lo SP	538°C	-73°C to 760°C (-100°F to 1400°F)
Alarm deviation*	Relay 1 AlmDevBand	3°C	1 to 56°C (1 to 100°F)
Inhibit alarm*	Relay 1 ALMInhibit	On	En or Dis
Type of event**	Relay 1 EventType	Time	Time or Temp
Event setpoint ***	Relay 1 Event SP	SPUpLim	SPLoLim to SPUpLim****

(repeat for relay outputs 2-4, if available)

Table of Settings

Relay	Туре	Action	Alarm/ Event Type	Setpoint	Alarm Deviation	Alarm Inhibit

^{*} appears only for alarm types

^{**} appears only for time or temperature event types

^{***} appears only for temperature event types

^{****} see enable page

^{*****} Normally de-energized and non-latching, normally energized and non-latching, normally de-energized and latching, normally energized and latching

Test

Menu Item	Display	Default	Range	Setting
Heater output	Test HeatOut	Off	On	
High limit relay	Test HiLimOut	Off	On	
Relay 1 output	Test Rly1 Out	Off	On	
Relay 2 output	Test Rly2 Out	Off	On	
Relay 3 output	Test Rly3 Out	Off	On	
Relay 4 output	Test Rly4 Out	Off	On	
High Limit Sensor	Test HL Temp	(sensor reading)		

Zone Cal

Menu Item	Display	Default	Range	Setting
Zone 1 actual	Zone Cal Zone1Act	38°C	-73°C to 760°C (-100°F to 1400°F)	
Zone1 displayed	Zone Cal Zone1Dis	38°C	-73°C to 760°C (-100°F to 1400°F)	
Zone 2 actual	Zone Cal Zone2Act	260°C	-73°C to 760°C (-100°F to 1400°F)	
Zone2 displayed	Zone Cal Zone2Dis	260°C	-73°C to 760°C (-100°F to 1400°F)	
Factory calibration*	Zone Cal FactCal	Ready	Ready or Done (push- key)	

^{*}only use when no calibration instrument is available

Sensor Cal

Menu Item	Display	Default	Range	Setting
Control Sensor Cal	Ctrl Sens 0F	-100°F	-100 to 1400°F	
	Ctrl Sens 1000F	1400°F	-100 to 1400°F	
	Ctrl Sens Done	Done	(read only)	
HiLim Sensor Cal	HL Sens 0F	-100°F	-100 to 1400°F	
	HL Sens 1000F	1400°F	-100 to 1400°F	
	HL Sens Done	Done	(read only)	
Retransmit Cal	RetCalLo XXXX *	0	0 to 4096**	
	RetCalHi XXXX *	4096	0 to 4096***	

^{*} note that the actual RetCalLo and RetCalHi values displayed are of no importance.
*** press ▲▼ keys until retransmission output = 1VDC
**** press ▲▼ keys until retransmission output = 5VDC

Enable Page

Menu Item	Display	Default	Range	Setting
Profiles	Enable Profile 1-8	Yes	Yes or No	
Autostart	Enable Auto Start	No	Yes or No **	
PID	Enable PID	Yes	Yes or No	
Control	Enable Control	No	Yes or No	
Communication	Enable Communication	No	Yes or No **	
Real Time Clock	Enable Clock	No	Yes or No **	
Relay outputs	Enable Relay 1-4	No	Yes or No **	
Test	Enable Test	No	Yes or No	
Zone Calibration	Enable Zone Cal	No	Yes or No	
Sensor Calibration	Enable Sensor Cal	No	Yes or No	
Setpoint lower limit	Enable SPLowerLim	20°C	-73°C to 759°C (-100°F to 1399°F)	
Setpoint upper limit	Enable SPUpperLim	260°C	-73°C to 760°C (-100°F to 1400°F)	
High limit overhead	Enable HiLimOH	5°C	3 to 11°C (5 to 20°F)*	
Password level 1	Enable Password 1	1	0 to 999	
Password level 2	Enable Password 2	2	0 to 999	
Manual/Timer mode setpoint change	Enable SPChange	Yes	Yes or No	
Analog Output Type	Enable Analog Type	Proc	Ctrl or Proc	

^{*} Maximum high limit setpoint (MaxHiLimSP) = SPUpperLim + HiLimOH Minimum high limit setpoint (MinHiLimSP) = SPLowerLim + HiLimOH

^{**} If the controller does not have this feature, "No" is the only option.

Technical Specifications

UL, cUL listed: UL file E136675

CE compliance to: • EMC Directive 89-366/EEC

European Standard EN55011/1991European Standard EN50082-2/1995

Power supply: • 100 to 240 VAC +10% -15%, 50-60Hz, 30VA Maximum

• 12 to 24 VAC/VDC +/-10%, DC to 60Hz, 30VA Maximum

Temperature: • Storage -20 to 60°C

Operating 0 to 50°C

Humidity: 90% or less, non-condensing

Sensor inputs: • Type J thermocouple -73°C to 760°C (-100°F to 1400°F)

• Input impedance 1M ohm or greater

Common mode noise rejection of 140db@60HzCommon mode input voltage of +/-12.0VDC

• Sample rate of at least 1 sample per second

• Stability of +/- 0.5°C per 5°C change in ambient temperature

 $\bullet~$ Repeatability of +/- 0.5 $^{\circ}\text{C},$ or +/-0.1% of sensed temperature (whichever is

greater)

Accuracy (@ 77°F +/-0.2 percent of span (+/-3°F)

Supply Voltage Influence of +/-0.5°C per 10% change in nominal line voltage

Temperature display:

1 degree resolution (C or F)

Accuracy after calibration of +/- 1°C, or +/-0.2% of sensed temperature

(whichever is greater) @25°C

• Four-digit seven-segment LCD, 0.43 inches high (11mm)

• Readout Stability (+/-1°F/10°F Change in ambient temperature)

Message display: • Two-line, 16 alpha-numeric 5x7 dot matrix characters per line

• 0.2 inches high (5mm)

Time base: +/- 4 seconds accuracy in 24 hours

Heat control output:

SSR Drive 24VDC nominal @ 70mA

Relay output: (plug-in

module)

Form A dry contact, rated 3 amps @ 24-264VAC

Retransmit output: 1 to 5 VDC into > 100K ohm load

Communications: (plug-in module)

RS232 Single drop, isolated
RS422 Multi-drop, isolated
RS485 Multi-drop, isolated

Remote inputs:

Dry-contact closure type with less than 250 ohm ON resistance

(plug-in module)

Front panel: NEMA 4X (with gasket)

Dimensions: 3.57H x 7.24W x 2.84D inches (91H x 184W x 72D mm)

Panel opening: 3.63H x 7.30W inches (92H x 185W mm)

MAINTENANCE

Do not attempt any service on this oven before opening the main power disconnect switch.

Checklist

- Keep equipment clean. Gradual dirt accumulation retards air flow. A dirty oven
 can result in unsatisfactory operation such as unbalanced temperature in the
 work chamber, reduced heating capacity, reduced production, overheated
 components, etc. Keep the walls, floor and ceiling of the oven work chamber
 free of dirt and dust. Floating dust or accumulated dirt may produce unsatisfactory work results. Keep all equipment accessible. Do not permit other
 materials to be stored or piled against it.
- <u>Protect controls against excessive heat.</u> This is particularly true of controls, motors or other equipment containing electronic components. Temperatures greater than 51.5EC (125EF) should be avoided.
- <u>Establish maintenance & checkup schedules</u>. Do this promptly and follow the schedules faithfully. Careful operation and maintenance will be more than paid for in continuous, safe and economical operation.
- <u>Maintain equipment in good repair.</u> Make repairs immediately. Delays may be costly in added expense for labor and materials and in prolonged shut down.
- <u>Practice safety.</u> Make it a prime policy to know what you are doing before you do
 it. Make CAUTION, PATIENCE, and GOOD JUDGEMENT the safety
 watchwords for the operation of your oven.
- <u>Lubrication</u>, Fan motor bearings are permanently lubricated. All door latches, hinges, door operating mechanisms, bearing or wear surfaces should be lubricated to ensure easy operation.

Tests

Tests should be performed carefully and regularly. The safety of personnel as well as the condition of equipment may depend upon the proper operation of any one of the functions of the temperature control. Test the control every 40 hours. Check that the heater LED is cycling on and off, indicating that the heater is working. Also check the high limit function to make sure it is working properly.

To test the high limit:

- 1. Go to Manual Mode and enter control setpoint value at least 20°F (11°C) lower than the current process temperature.
- 2. Press the Menu key and lower the high limit setpoint to a value just below the current process temperature.
- 3. The high limit alarm indicator should flash and a high limit alarm message should be displayed. Verify that the heater relay has been disabled by checking that 2LED on the control panel circuit board is not lit.
- 4. Return the control setpoint and high limit setpoint values to their original values.
- 5. Press the Reset key.

Nitrogen Supply

Turn the three-way valve on the control panel to the OFF position and screw the adjusting knob clockwise on both PURGE and MAINTAIN flowmeter all the way to the OFF position. Then open the tank valve on the nitrogen supply and set the pressure regulator to about 40 PSI. Check the nitrogen plumbing for leaks

WARNING:

Disconnect the main power switch or power cord before attempting any repair or adjustment.

with a soapy water solution. As nitrogen gas is odorless, all leaks should be stopped to prevent the possibility of suffocation in a small work areas. Gas from a nitrogen leak might displace much of the oxygen in the atmosphere.

Replacement

Parts

To order or return parts, contact the Service Products Division at Despatch. The Service Products features our Response Center for customer service. When returning parts, a Despatch representative will provide you with an MRA (Material Return Authorization) number. The MRA number must be attached to the returned part for identification. When you are ordering parts, be sure to give the model number, serial number and the part number. This will expedite the process of obtaining a replacement part.

When you have a **service** need, contact the **Response Center** at **1-800-473-7373**: FAX 612-781-5353.

WARNING:

Disconnect the main power switch or power cord before attempting any repair or adjustment.

Protocol Plus™ Instrument

(Tools needed: 1/4" socket driver with #1 bit, #2 Phillips screwdriver)

- 1. Disconnect the power.
- 2. Remove the screws from the control panel and slide it forward.
- 3. Unplug all terminals on the rear of the control, noting the proper connections.
- 4. Remove the four retaining clips for the control.
- 5. Remove the control.
- 6. Insert the new control into the panel.
- 7. Fasten the four retaining clips.
- 8. Re-plug all terminals.
- 9. Fasten the control panel.

Heater Unit

(Tools needed: 3/8" wrench, square recess driver with #1 bit for LND1-42, one-quarter inch socket set for LND2-11 and LND2-24)

- 1. Remove the chamber floor plate.
 - a. Remove the screws from the floor plate.
 - b. Lift the floor plate out of the oven.
- 2. Disconnect the heater leads from heater element with wrench. Note which wires go on which terminals.
- 3. Unscrew the screws holding the heater frame to the oven body.
- 4. Remove the heater and discard.
- 5. Screw down the new heater frame.
- 6. Attach the heater leads to appropriate terminals.
- 7. Replace the interior floor and screws.

Fan Motor

(Tools needed: Screwdriver, 5/32 inch Allen wrench, one quarter (1/4) inch socket set)

- 1. Remove the chamber floor plate.
 - a. Remove the screws from the floor plate.
 - b. Lift the floor plate out of the oven.
- 2. Remove the left side wall.
- 3. Remove the fan inlet plate.
- 4. Loosen the set screws (2) on fan wheel in middle of oven bottom. You can reach the fan wheel by going through the heater or by disconnecting and removing the heater. Refer to the Heater Unit instructions.
- 5. Place the oven on its back (LND1-42 only).
- 6. Remove the bottom plate on the LND1-42, or remove the control panel on the LND2-11 and 2-24. This will reveal the fan motor.
- 7. Remove the fan motor.
 - a. Unscrew the screws (4) holding motor mounts to the body.
 - b. Disconnect the motor leads from the terminal strip on the power panel.
 - c. Lift the fan motor from the oven body.

After the fan wheel has run at temperature for awhile, it will stick to the shaft. Some force may be required to separate the fan wheel from the fan motor shaft.

- 8. Take the motor mounts off old motor.
- 9. Put the motor mounts onto new motor.
- 10. Replace the fan motor.
 - a. Insert the shaft into shaft collar. Put the fan wheel onto shaft from inside oven.
 - b. Reattach the motor mounts to the oven body, making sure grommets are in place.

- c. Reattach the motor lead wires to the terminal strip.
- 11. Replace the oven bottom (LND1-42) or control panel front (LND2-11 or 2-24).
- 12. Turn the oven right side up.
- 13. Adjust the fan wheel for 3/16 inch clearance between the wheel and the inlet ring.
- 14. Tighten the set screws on the fan wheel.
- 15. Check that the set screws hit the flats machined into the motor shaft.
- 16. Replace the chamber floor plate.

TROUBLESHOOTING

Equipment which operates for long periods of time may develop problems. Below are possible problems and suggested solutions. If you have a problem not listed and do not know what to do, contact Despatch Industries at our toll free Help Line 800-473-7373.

Difficulty	Probable Cause	Suggested Remedy
Failure to heat	No power	Check power source and/or oven and wall fuses.
	Burned out heater	Replace heater (see warranty.)
	Protocol™ malfunction	Replace controller.
	Loose wire connections	Disconnect power and check connections behind control panel.
	Purge timer not timed out or failed	Wait 3 minutes. If still not heating, replace purge timer.
	Heater relay failure	Replace circuit board.
	Airflow switch not closed	Check exhaust system for proper operation and ventilation. If airflow switch still does not close, replace switch.
	Door switch failure (if installed)	Replace door switch.
Slow heat up	Improperly loaded	Reduce load or redistribute load in chamber.
	Low line voltage	Supply sufficient power and proper connections. Check for circuit overload.
	Heating element(s) are burned out	Replace burned out element (see warranty statement.)
	240 volt oven is connected to a 208V line (LND1-42 only)	Raise line voltage to a 240 volt line or modify oven for 208V operation (consult factory).
	Fan motor failure	Replace fan motor.
Frequent heater element out	Harmful fumes generated by load	Increase vent opening or discontinue process.
	Spillage or splattering of material on heater elements	Disconnect power and clean oven chamber and elements.
	Overheating oven	Check the Hi-limit.

Difficulty	Probable Cause	Suggested Remedy	
Erratic temp. or inaccurate temp.	Protocol™ malfunction	Replace Protocol™.	
	Improper tuning parameters	Check tuning parameters.	
	Protocol™ miscalibration	Recalibrate Protocol™ (see section on Calibration mode.)	
	Hi-limit setting	Hi-limit should be 10-25EC higher than setpoint.	
	Improper offset	Check zone calibration.	
Excess surface or door temp.	Door seal deterioration	Replace door seal.	
Improper airflow	Fan motor failure	Replace fan motor.	
	Fan wheel seated too low on fan shaft	Adjust fan wheel for 3/16" clearance between wheel and inlet ring.	
	Unbalanced fan wheel	Replace fan wheel.	
Excessive vibration	Dirty fan wheel	Clean fan.	
	Unbalanced fan wheel	Replace fan wheel.	
Oven will not control at	Hi-limit set too low	Set the Hi-limit higher	
setpoint	Protocol malfunction	Replace control.	
	SSR malfunction	Replace SSR and/or check control output voltage.	
		Open exhaust air vent. Unit will not control below	
	Air friction of recirculation fan	minimum operating temperature with vent closed.	
Heater does not	Protocol malfunction	Replace Protocol	
shut down until	000 " "		
temp. reaches the Hi-limit	SSR malfunction	Replace SSR	
setting			

NOTE: The circuit board mounted on the control panel has three status LED indicators to help troubleshoot if the oven is not heating (LND1-42 only).

- A. If LED 1 is not lit, check 2F and 3F (control fuses), or power switch.
- B. If LED 1 and LED 3 are lit but not LED 2, check high limit (and optional door switch, if installed).
- C. If all three LEDs are lit, check 1F and 4F (heater fuses), SSR, heater, and heater relays.

APPENDIX

Temperature Scale Conversion (C/F)

The Protocol Plus controller can be operated in either C or F. The default setting for the controller is C. Changing from one to the other is as follows:

- 1. Go into the **Setup Mode** on the controller.
- 2. Press the **Select** Key until **Setup** is displayed.
- 3. Press the **Page** key and **Security** will be displayed.
- 4. Press the **Menu** Key and **Password** will be displayed. Use the **arrow** keys to enter the proper password. The default password is 2 for level two.
- 5. Once the proper password is displayed, press the **Page** key until **PID** is displayed.
- 6. Press the **Menu** key and **Temp Unit** along with C or F will be displayed. Use the arrow keys to change the setting.
- 7. Once the proper setting is displayed, press and hold the **Page** key for approximately three seconds to exit the **Setup Mode**.

Optional MRC5000 Recorder Setup

The temperature is retransmitted to the Recorder from the Controller. Set up the Recorder as follows:

- 1. Make sure that jumper **JU1** is setup for the **5 VDC** setting (see MRC Manual).
- 2. Move the **Mode** switch to the **PROG/TEST/CAL** position, and **Prog** will be displayed.
- 3. Press the **down arrow** key twice and **Inps** will be displayed. Make sure the settings are per the table below.
- 4. Once all the settings have been changed, move the **Mode** switch to the **RUN** position. The display on both the Recorder and controller should read the same.

Parameter Code	Degrees C	Degrees F
Inps	17	17
Icor	0	0
diSP	On	On
dPOS	0	0
EUU **	400	752
EUL **	0	32
ChUP	400	800 *
ChLO	0	0
DFF	1	1

* Note: The 0 - 400 chart paper must be changed to the 0 - 800 chart paper. Depending on the equipment, 0 - 600 paper can be used if the maximum temperature is 500 degrees F.

** Note: These values must mach the setting sRetOutLo and RetOutHi on the Control page on the Protocol Plus controller (example: RetOutLo is 32, then EUL must read 32).